# Deep Freezer Gudero+ Series

Users Manual, Ver. 4.0

30/03/17

**Failure to read, understand and follow the instructions in this manual may result in damage to the unit, injury to operating personnel and poor equipment performance.**

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These operating instructions are protected by copyright. Rights resulting thereof, particularly reprint, photomechanical or digital post processing or reproduction, even in part, are only allowed with the written consent of ilShinbiobase.

This regulation does not apply to reproductions for in-plant use.

## **Trademarks**

All trademarks mentioned in the operating instructions are the exclusive property of the respective.

ilShinbiobase Co. Ltd. provides this document to its customers with a product purchase to use in the product operation. This document is copyright protected and any reproduction of the whole or any part of this document is strictly prohibited, except with the written authorization of ilShinbiobase Co. Ltd.

The contents of this document are subject to change without notice. All technical information in this document is for reference purposes only. System configurations and specifications in this document supersede all previous information received by the purchaser.

ilShinbiobase Co. Ltd. makes no representations that this document is complete, accurate or error-free and assumes no responsibility and will not be liable for any errors, omissions, damage or loss that might result from any use of this document, even if the information in the document is followed properly.

This document is not part of any sales contract between ilShinbiobase Co. Ltd. and a purchaser. This document shall in no way govern or modify any Terms and Conditions of Sale, which Terms and Conditions of Sale shall govern all conflicting information between the two documents.

**© ilShinbiobase Co. Ltd. All rights reserved.**

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**(Read and understand the manual before using this machine)**

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**(While machine works, the printed sign can be detached with poor adhesion, then attach with the same one)**

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**Only authorized personnel may service this equipment**

**Before Installation**

## -Transportation-

If the unit is tilted excessively, the compressor lubricant may flow from the compressor (or oil separator) to the line system. Insufficient filling level may damage the compressor when the refrigeration system is started. Keep vertical position while transporting the unit. During transportation, the unit must not tilt more than 10° .

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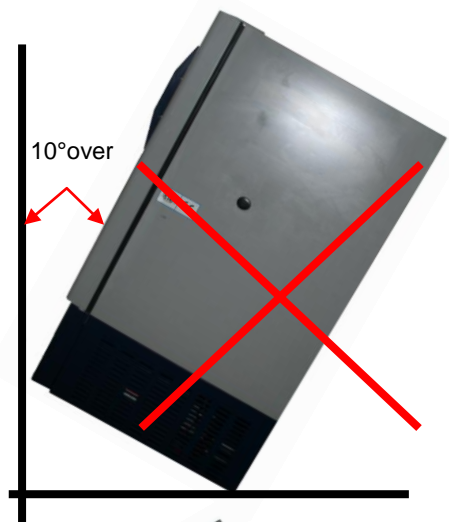
Max. tilt angle transportation (Enable position)

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
### **1. Recommended room condition (권장 장소)**

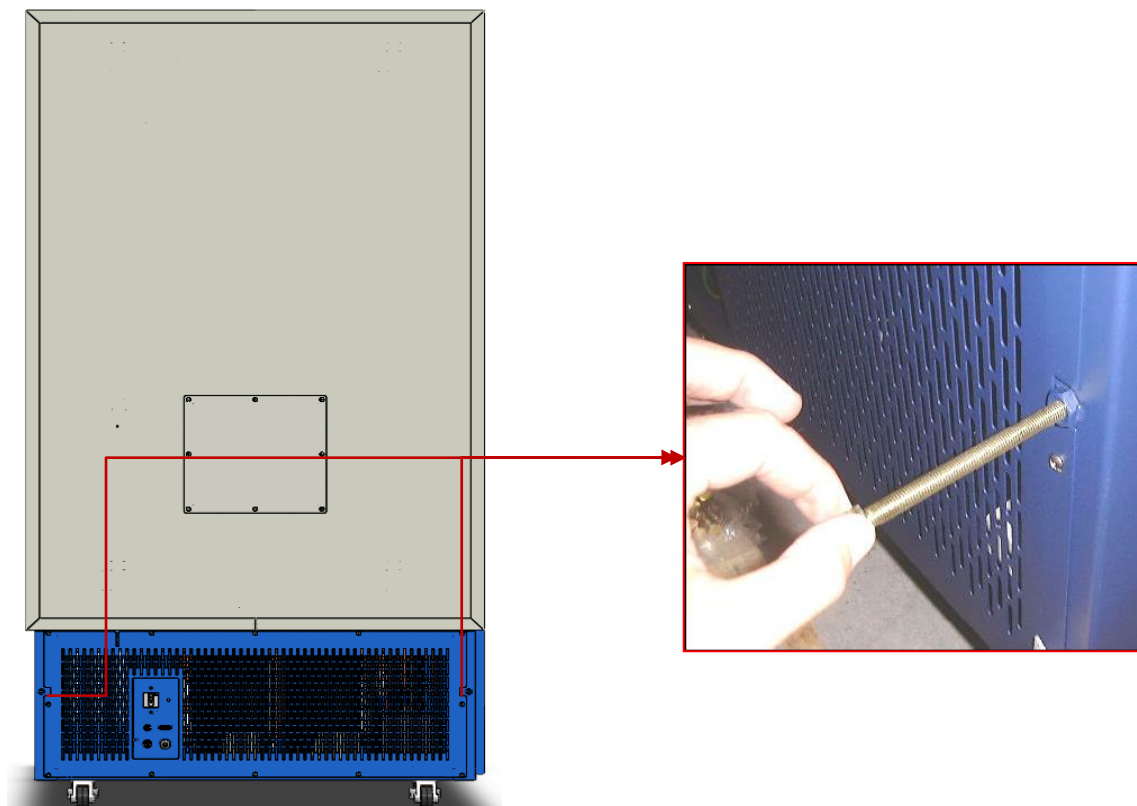
- a. Recommended room temperature: between 5℃ and 28℃. If the room temperature is over 30℃, please ensure that suitable air-conditioning is provided.
  - b. Make sure the floor is level and even.
  - c. Make sure the condenser grille is not blocked. (Adequate air circulation is required)
  - d. Install the two spacer pins on the back of the freezer to allow heat dispersion behind the freezer.
- a. 권장 대기 온도: 5℃ ~ 28℃ 만약 실온이 30℃보다 높으면, 적절한 냉난방시설이 갖추어 지기를 권장합니다.
- b. 바닥은 수평이고, 평평한 곳에 설치해 주십시오.
- c. 장비는 충분한 공기 순환이 필요합니다.
- d. Freezer 뒤로 열이 배출되도록 freezer뒤에 두 개의 spacer pin을 설치해 주십시오.

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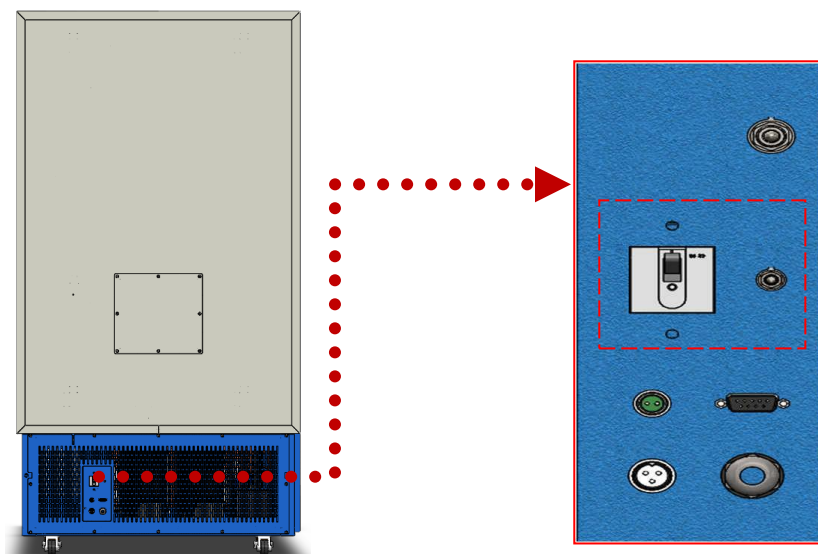**We wish you a quick and trouble-free installation!**

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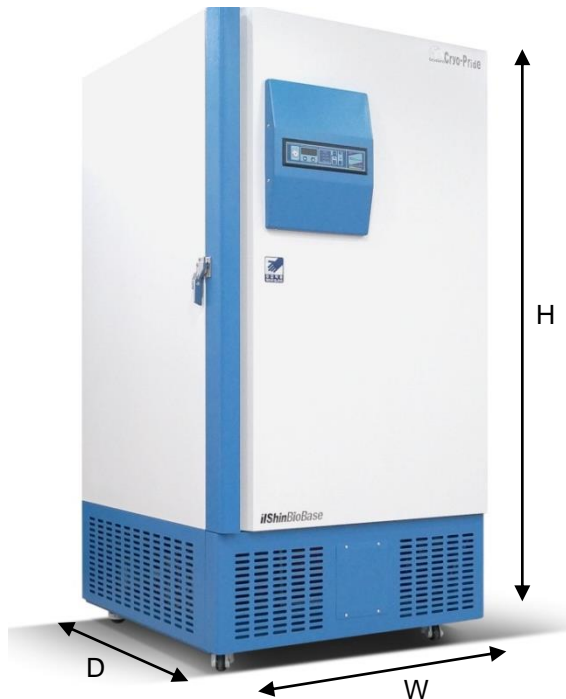
## 1. Dimension of The Freezer Model (모델별 규격)

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
Model No.	External Dimensions	Internal Dimensions	Capacity(L)	Weight(kg)
	WxDxH	WxDxH		
DF3510	920x805x1660	600x480x970	276	256
<b>DF8510</b>				
DF3514	920x805x1980	600x480x1290	369	279
<b>DF8514</b>				
DF3517	920x955x1980	600x630x1290	484	307
<b>DF8517</b>				
DF3520	1070x955x1980	750x630x1290	605	340
<b>DF8520</b>				
DF3524	1227x967x1980	890x635x1290	730	378
<b>DF8524</b>				
DF3530	1227x985x1980	890x725x1290	832	410
<b>DF8530</b>				

The operating temperature of Model DF35\*\* is  $-10^{\circ}\text{C}$  to  $-35^{\circ}\text{C}$ , and for the Model DF85\*\* this is  $-60^{\circ}\text{C}$  to  $-90^{\circ}\text{C}$ . (normal temperature for usages is  $-20^{\circ}\text{C}$  to  $-30^{\circ}\text{C}$ ,  $-60^{\circ}\text{C}$  to  $-80^{\circ}\text{C}$ ). In addition, if the last two numbers(\*\*) of the model are the same, the same operating temperature is applicable.

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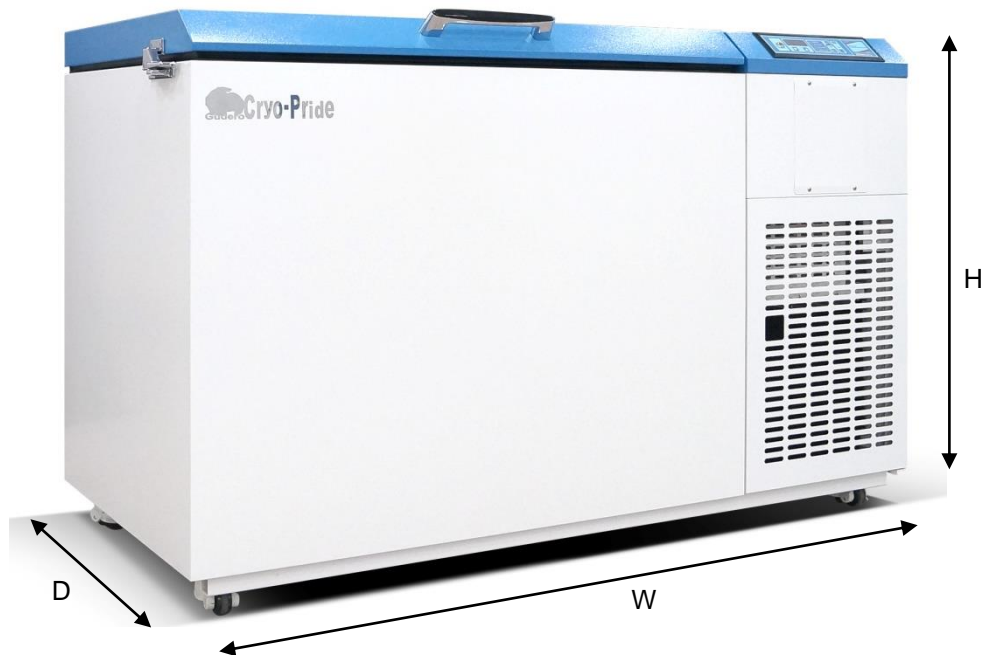The operating temperature of Model DF45\*\* is -10℃ to -45℃, and for the Model DF\*\* 90 -40℃ to -90℃ (normal temperature for usages is -20℃ ~ -40℃, -60℃ to -85℃). In addition, if the last two numbers(\*\*)of the models are the same, the same operating temperature is applicable.

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
## 2. Specification (기술 사양)
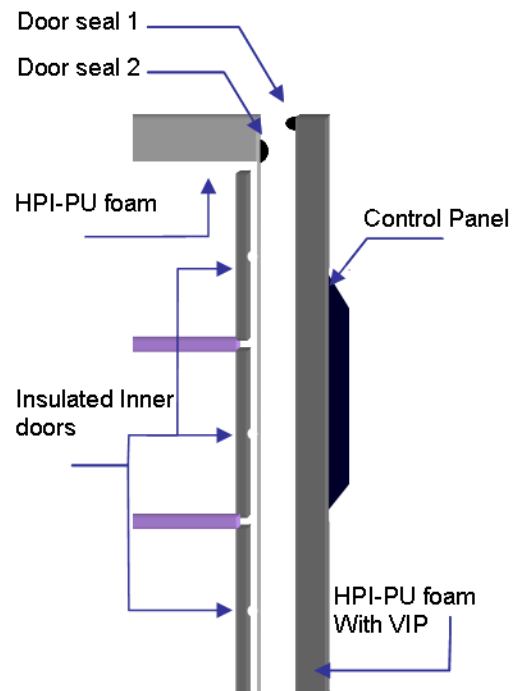
### 2.1 Deep Freezers Characteristics (초저온 냉동고 특징)

- I. Implementation of durable hermetic compressors
- II. ECO mode considering power consumption with better efficiency
- III. Efficient evaporator with minimum contact for fast cooling down of chamber
- IV. 130mm polyurethane foamed insulation with VIP® insulation
- V. Heavy duty casters with built in adjustable stands(unique)
- VI. Cabinet with round corners for easy cleaning
- VII. 2 built-in access ports
- VIII. User-friendly control panel to eye level
- IX. USB data download port and dust-free condenser design
- X. Pressure equalization port allows easy access to cabinet after door closing
- XI. Quiet running fans and compressors create very low noise level (50~65 dB(A))
- XII. Heated door seals with hot gas prevent ice build-up

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
#### Theory Behind the Perfect Fit System

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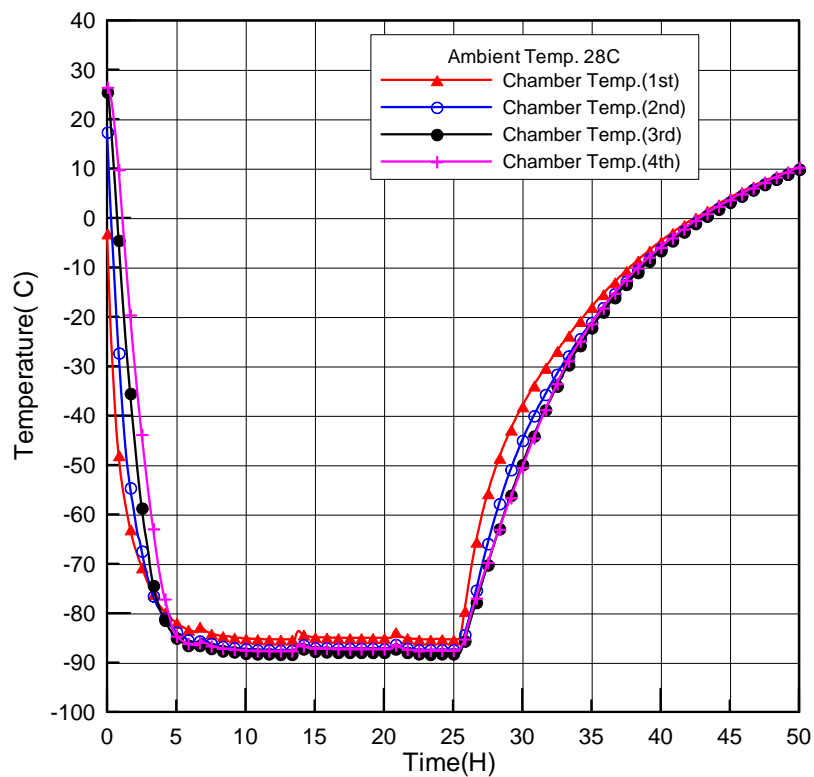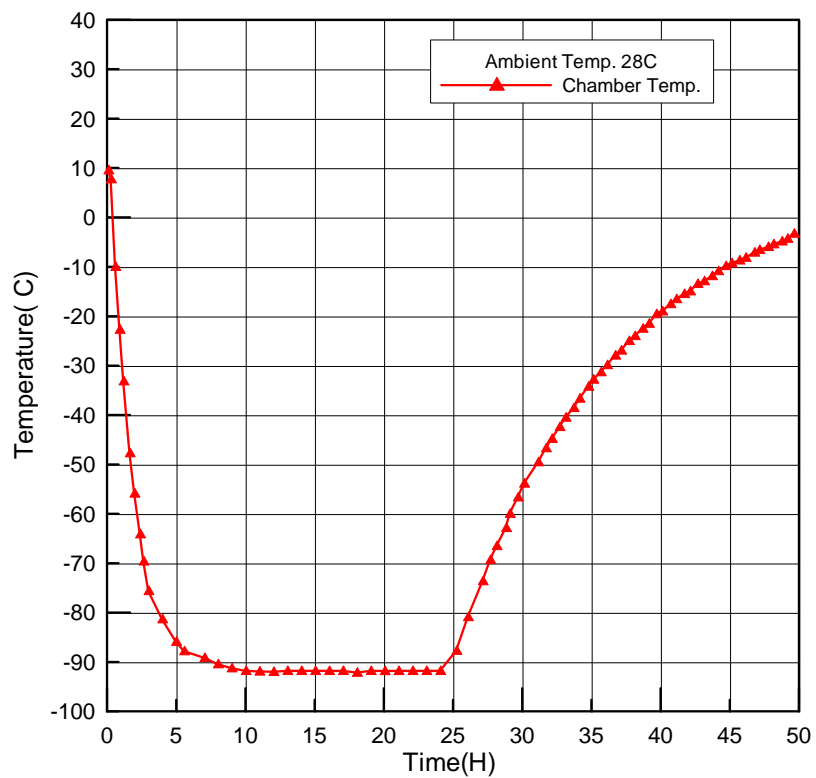
## Installation (설치)

### 3.1 Installation Place (설치 장소)

1. Avoid direct sunlight
2. Adequate air circulation
3. Not subject to excessive dust
4. Not subject to mechanical vibration or shock
5. Ambient temperature 5~33℃, optimum temp. 18~ 25℃
6. Operating Height : Below 2,000m
7. Operating moisture: Below 80% RH, if the max. operating temperature is 33℃, it should be below 57% RH
8. Supply voltage fluctuation : Lower than 10%

\* Caution: If you install the unit other places than mentioned, the unit may not operate properly.

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### 3.2 How to Install (설치 방법)

1. Keep the distance of 15cm or longer between the wall and the backside of the unit.
2. Do not install the unit in an uneven position.
3. Use castor fixing nuts on the slippery floor.
4. Condenser grille should not be obstructed

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1. To ensure secure connection, connect the power plug firmly to the power socket.
2. Don't connect the plug of the unit together with plugs of other units in the same socket.
3. Make sure that the power supplying cord is not folded and pressed.
4. If you use an extension wire, it must have capacity above 250V, 15A and not longer than 3m.
5. This unit is exclusively designed for 220(230)V, 60(50)Hz, 30Amp. Please check the operating voltage before use. For 110V, use transformer(above 3kW).

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### 4.2 How to Earth (접지 방법)

1. For security reasons, the unit should be earthed.
2. If the power socket is fitted with earth rod (230V), please check the earth terminal before use.
3. If the power socket is not fitted with earth rod, connect the earth wire to the unit. Then, connect it to the copper plate. Finally, bury it in the damped ground(25cm under the ground or deeper).

\* Warning: Do not earth to gas pipe, telephone line or lightning rod.

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The rating Amp. of JABA Series circuit breaker must be equal or higher than 30. Please check before installation.

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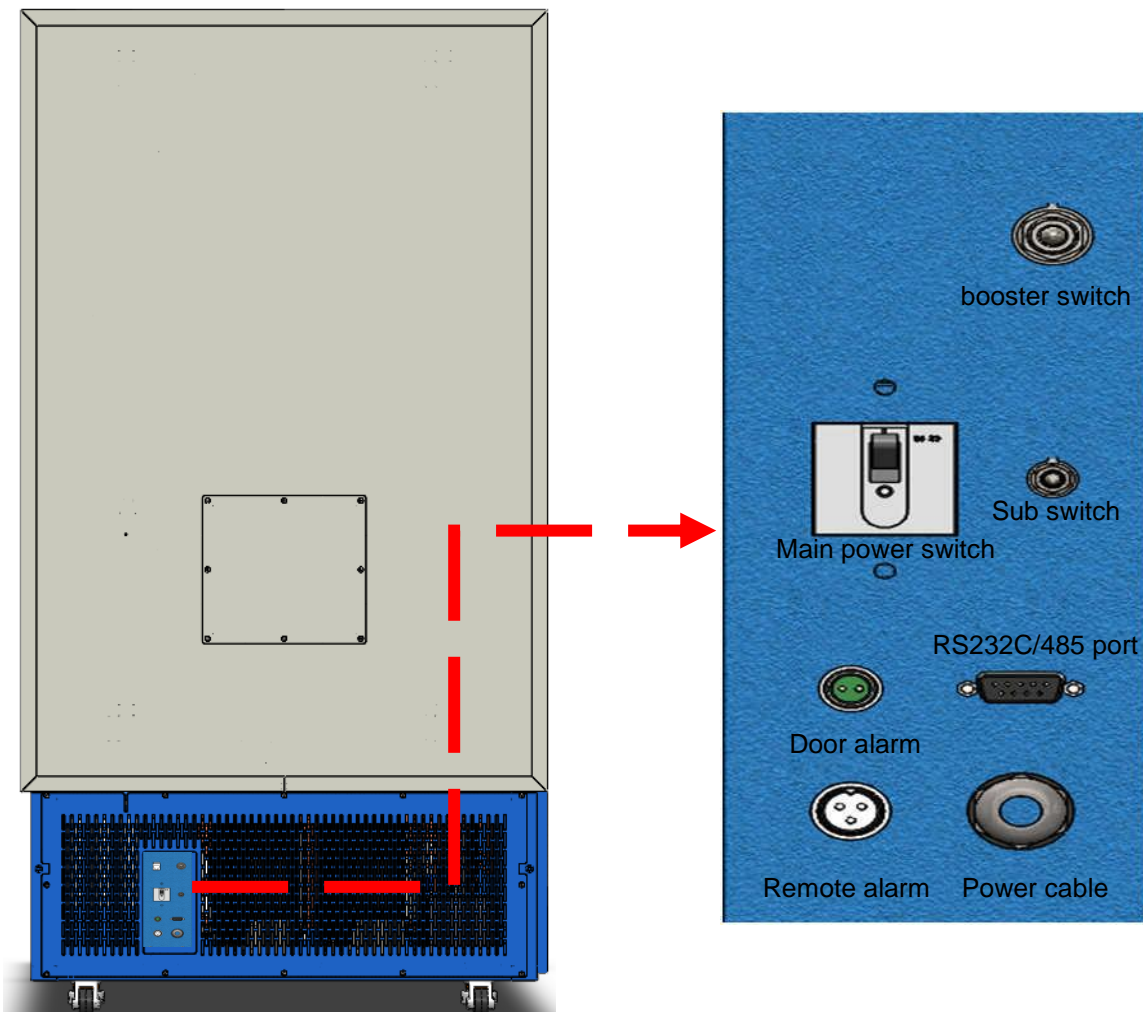
(Main power switch is located on rear side of the machine. User must turn on “Main power switch” first and turn on “Sub switch”. When user needs to turn off the machine, turn off “Sub switch” first and turn off “Main power switch”. When power is connected, touch screen will be activated.

**\*Notice : Circuit is not protected, when blackout happens, during operation by Main power switch**

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#### 4.4 Door Alarm (도어 알람)

The door alarm is used in combination with the CO<sub>2</sub> backup system. It is safe to make sure that the operator will not get liquid CO<sub>2</sub>/LN<sub>2</sub> in the face in case of opening the door of the freezer. Always connect the CO<sub>2</sub>/LN<sub>2</sub> backup system to the door alarm on the back of the freezer.

For safety, measure resistance on the contacts with Ohmmeter

1. Measure contacts of door alarm and resistance shows 0 Ohm. Door is closed.
2. Open the chamber door.
3. Measure contacts of door alarm again and resistance shows limitless. Door is open.

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## 4.5 Remote(External) Alarm (리모트(외부) 알람)

The remote alarm is a voltage free dry contact which will alarm in case of,

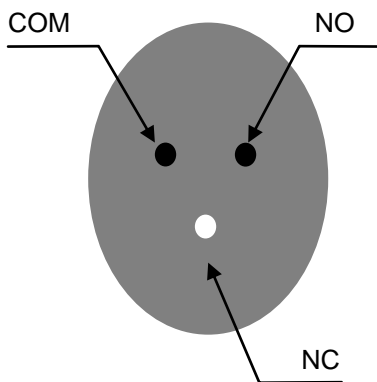
1. High temperature alarm
2. Low temperature alarm
3. Power failure
4. Clean filter alarm

Under normal condition, relay is always closed which means COM and NC (Normally Closed ) are connected. In case of alarm, relay is open and COM and NO (Normally Open) are connected.

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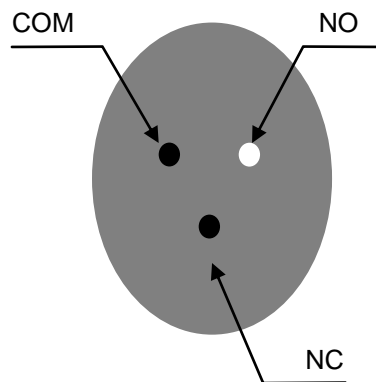
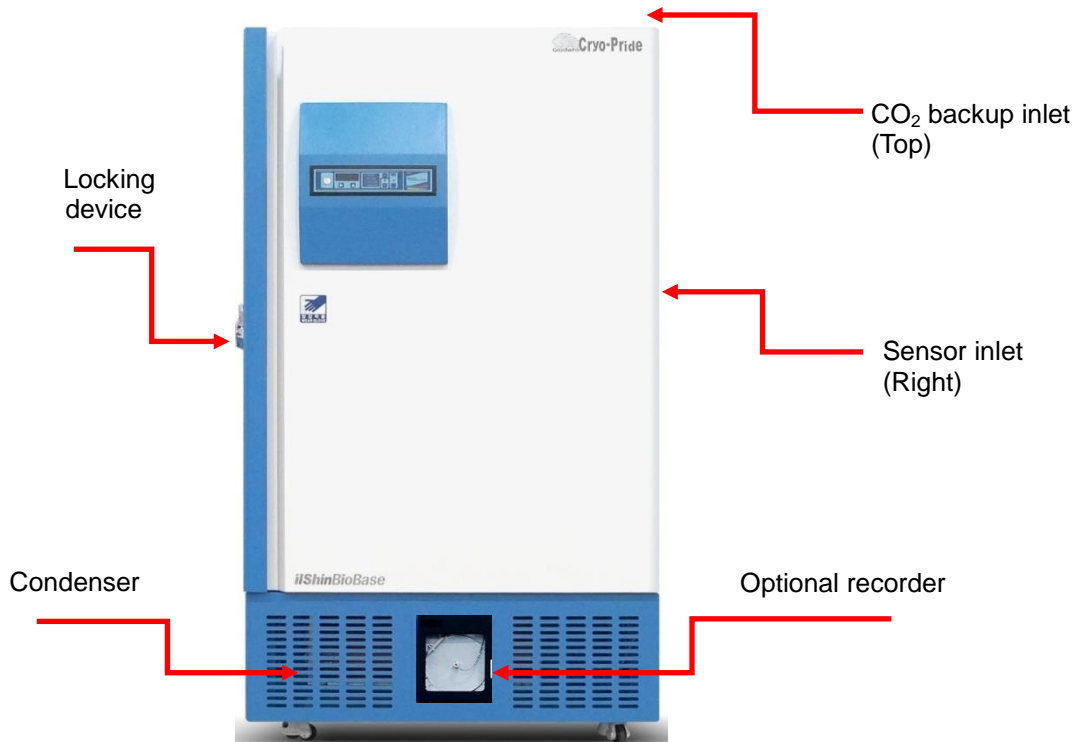
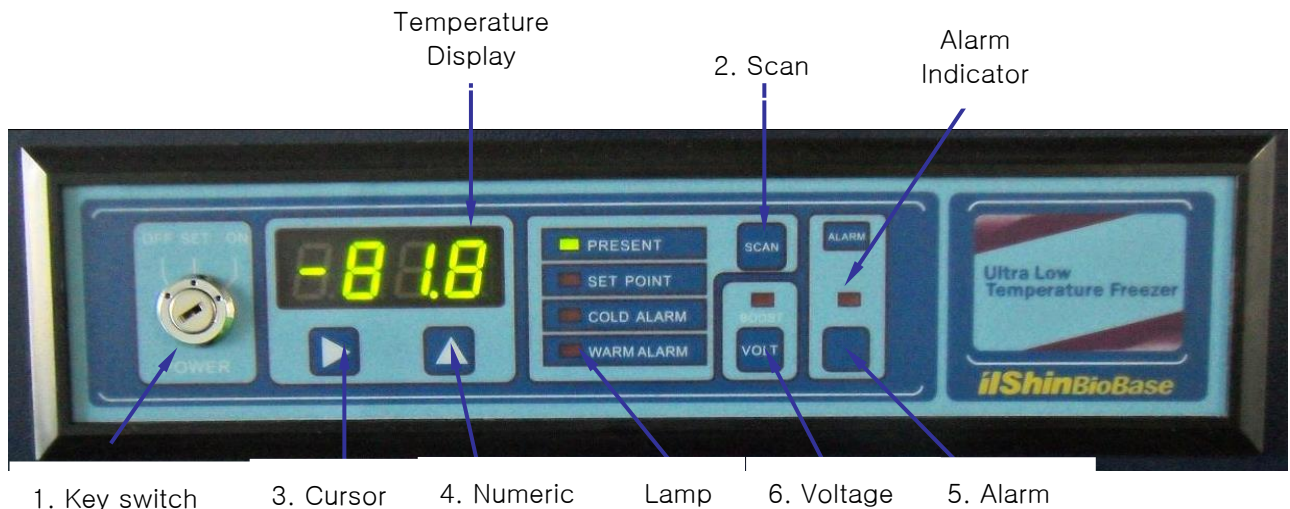
NO.	BUTTON	CONTROLS
1.	KEY SWITCH (키 스위치)	<ul style="list-style-type: none"> <li>- During OFF, all functions are stopped</li> <li>- During SET, all functions are working, and it is possible to adjust the temperature</li> <li>- During ON, the unit is operating as planned, and it is not possible to adjust the temperature.</li> </ul> <p>* Function blocking system : If you take out the key during ON, all functions will be blocked (Please keep the key well, for the planning of functions)</p> <ul style="list-style-type: none"> <li>- OFF상태에서는 모든 작동이 정지합니다.</li> <li>- SET상태에서는 모든 기능이 작동되며 온도수정이 가능합니다.</li> <li>- ON상태에서는 설정된 대로 정상 작동하며 온도 수정은 할 수 없습니다.</li> </ul> <p>* 기능잠금장치 : KEY를 ON상태에서 빼시면 기능 잠금상태로 됩니다. (다음 작동을 위해서 키를 잘 보관하십시오.)</p>
2.	SCAN (스 켄)	<ul style="list-style-type: none"> <li>- This button is to select each temperature mode (Only in the case that the KEY SWITCH is set to SET)</li> </ul> <p>- 각 온도 모드를 선택 및 설정하는 사용합니다. (KEY SWITCH를 SET로 하여 각 온도를 설정하고자 할 때 사용합니다.)</p>

--	--	--

NO.	BUTTON	CONTROLS
3.	CURSOR (위치변환)	<ul style="list-style-type: none"> <li>- This button is to select a position of numerical value</li> <li>- 변경하고자 하는 수치의 위치를 선택할 때 사용합니다.</li> </ul>
4.	NUMERIC (숫자변환)	<ul style="list-style-type: none"> <li>- This button is to change a numerical value</li> <li>- 수치를 변경할 때 사용합니다.</li> </ul>
5.	ALARM (알 램)	<ul style="list-style-type: none"> <li>- If the inside temperature of the unit exceeds the alarming temperature range, an error message will come on the temperature indicator. The lamp will be on and the alarm signal will sound. (Push the alarm button (mute), then, the alarm signal does not sound)</li> <li>- 고내온도가 상하한 경보온도를 벗어날 때 온도표시기에 에러 메시지가 표시되며,ALARM 램프가 점등, 경보음이 울립니다. (ALARM버튼을 누르면 경보음은 정지합니다.)</li> </ul>
6.	VOLTAGE (전 압)	<ul style="list-style-type: none"> <li>- This button indicates the voltage (Push the button, then the voltage will be indicated on the Temperature indicator</li> <li>- 현 전압을 확인하고자 할 때 사용합니다.(VOLT 버튼을 누르면 온도 표시기에 현 제품을 인입되는 전압이 표시됩니다.)</li> </ul>

## 5.4 How to Operate (작동 순서)

1. Turn the KEY SWITCH to the right and set it to SET.
2. All functions will be operational. The PRESENT lamp will be on and the inside temperature of the Chamber will be shown.
3. Push the SCAN button once to check the temperature. The SET POINT lamp will be on. (To adjust SET POINT, use the CURSOR button and NUMERIC button)
4. Push the SCAN button once more to check minimum alarm temperature. Then, the COLD ALARM lamp will be on. (To adjust the COLD ALARM temperature, use the CURSOR button and NUMERIC button) The COLD ALARM temperature should be set lower than the SET POINT.  
 \* Temperature : e.g. In the case that the SET POINT temperature is  $-70^{\circ}\text{C}$  , the COLD ALARM temperature is  $-75^{\circ}\text{C}$  .
5. Push the SCAN button once more to check the maximum alarm temperature. Then, the WARM lamp will be on. (To adjust the WARM ALARM temperature, use the CURSOR button and NUMERIC button)  
 \* The WARM ALARM temperature should be set higher than the SET POINT temperature:  
 e.g. In the case that the SET POINT temperature is  $-70^{\circ}\text{C}$  , the WARM ALARM temperature is  $-65^{\circ}\text{C}$  .
6. Push the SCAN button once more to turn on the PRESENT lamp showing the present inside temperature of the deep freezer.

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7. Push the SCAN button once more to turn on the PRESENT lamp showing the present inside temperature of the deep freezer.
8. Now, everything is set. Turn the KEY SWITCH to the right once more and set it to ON. Take out the KEY by pulling it to the front. Please keep the KEY carefully to change the temperature in the future. (As long as the key is taken out, it is impossible to set the temperature. It prevents the unit from changing the temperature by the third party or by mistake. In this way, the test material and the unit can be protected).
9. To check the voltage, push the VOLT button, then, the current incoming voltage will be shown on the temperature indicator.
10. **The voltage Booster is equipped :**
  - i. If the input voltage is lower than 205V, the lamp of the Booster will be on and the voltage boosted. **e.g. 205V->225V (15V boosted)**
  - ii. If the input voltage is lower than 180V, the operation of the unit will be stop to protect the parts of the unit. (The functions such as temperature indication and alarm, are still operating)

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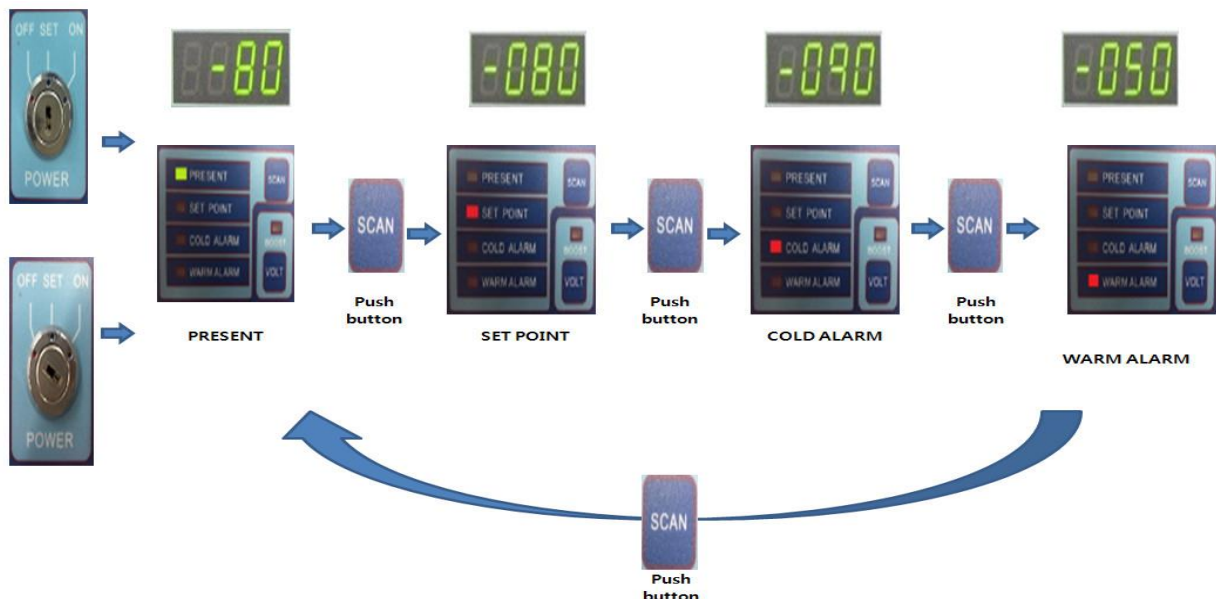
### 6.2 Fireboard Storage Boxes and Dividers (박스과 칸막이)

#### External dimensions:

External dimensions of 2 inch boxes (model PP0060): 133x133x49 mm (DxWxH)

External dimensions of 3 inch boxes (model PP0061): 133x133x77 mm (DxWxH)

#### Internal Dimensions:

Internal dimensions of 2 inch boxes (model PP0060): 124x124x45 mm (DxWxH)

Internal dimensions of 3 inch boxes (model PP0061): 124x124x70 mm (DxWxH)

#### **Possible dividers:**

Dividers are also made of plasticized card box. Height of all dividers is 35 mm

#### Model: Description:

PP1049 Storage box divider 7 x 7

PP1064 Storage box divider 8 x 8

PP1081 Storage box divider 9 x 9

PP1100 Storage box divider 10 x 10

Thick plasticized card box material

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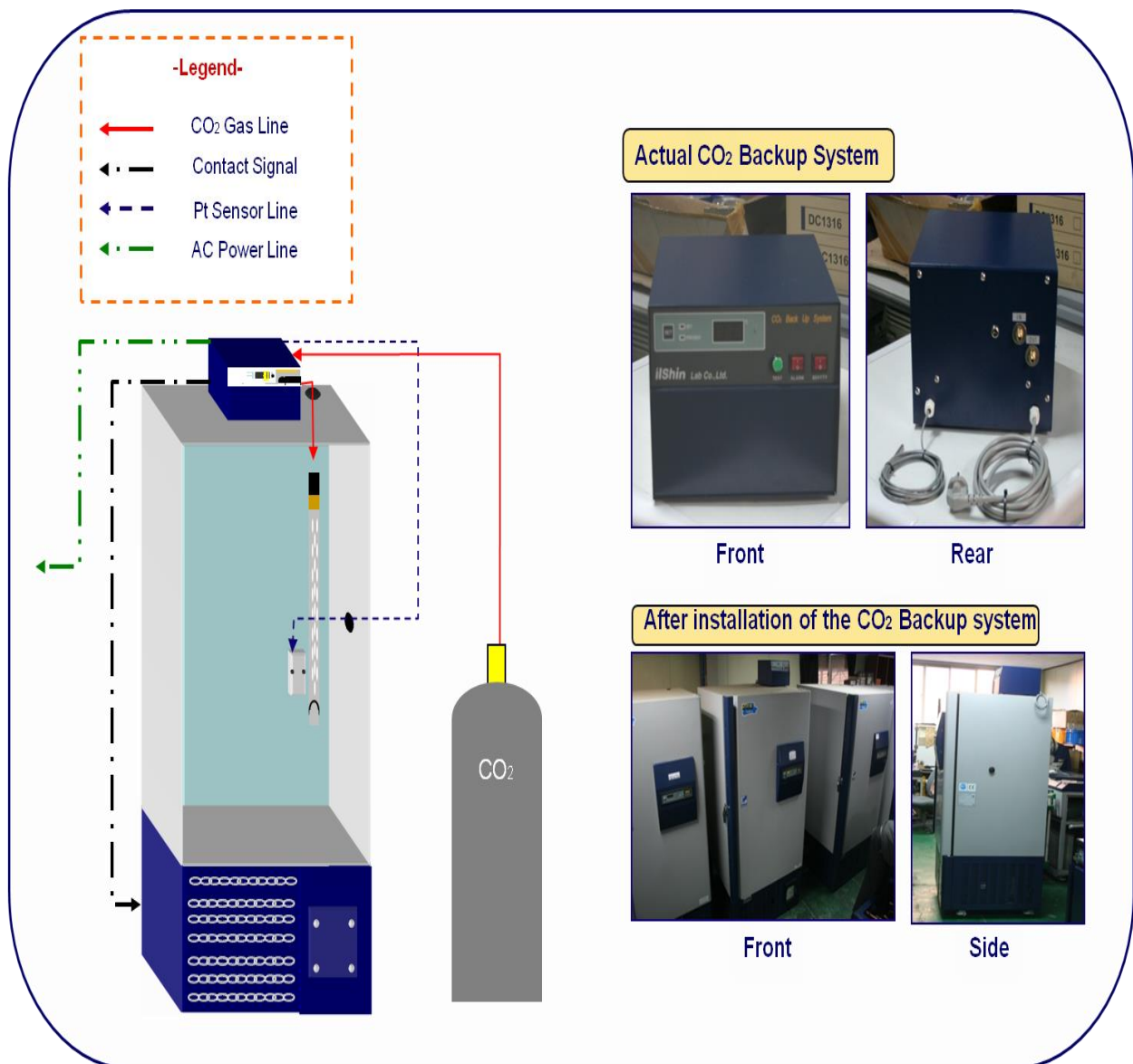
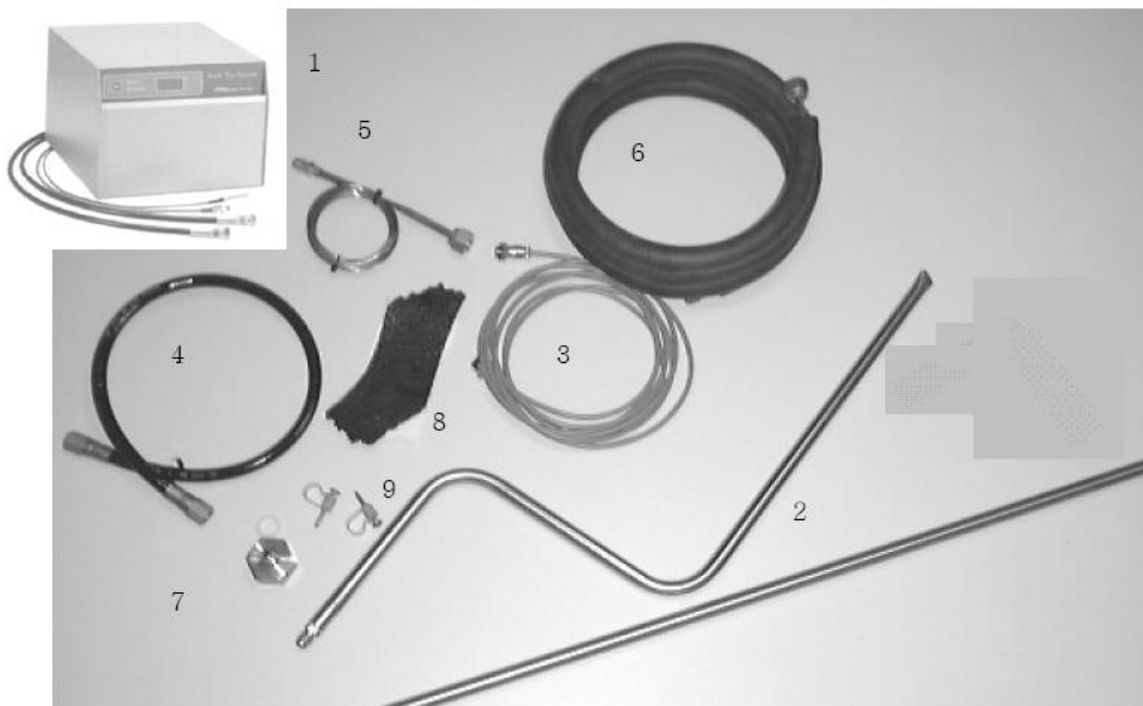
This backup system is used in case that the temperature goes up due to failure of power or technical problem. The temperature set to BACK UP SYSTEM will be observed. As a result, the damage of the sample in the unit is minimized.

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#### 6.3.1 Index of the parts (부분별 명칭)

< CO<sub>2</sub> backup system structure >

1. CO<sub>2</sub> backup system
2. Injection tube with hoses
3. Connection cable for door switch
4. Black flexible hose
5. Capillary tube
6. Copper insulated hose
7. Connection for CO<sub>2</sub> cylinder (including Teflon ring for better fitting)
8. Roll kit for sealing up
9. Connection screws (2x)

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1. Connect a temperature resistant tube between the opening of CO<sub>2</sub> cylinder and the ④ IN of the rear side.
2. Insert the injection rod of CO<sub>2</sub> into the inlet on top of the unit, and connect the rod to the ⑤ OUT of the rear side of Control System with a tube.
3. Insert the Sensor Line into the holes at the back and right side.
4. Open the valve of CO<sub>2</sub> cylinder and check by soap water whether there is a leakage around connection part.
5. Connect the power supply. (If the inside temperature of the deep freezer is higher than the set temperature, CO<sub>2</sub> gas will be injected. Please connect the power supply under the set temperature.)
6. Switch on the BATTERY SWITCH
  - \* N.B.: The BATTERY SWITCH should be OFF during moving, storage, or POWER OFF. (When the BATTERY SWITCH is on for a long time with power OFF, the battery will be discharged and should be replaced.)
7. Switch the ALARM SWITCH on, to activate the alarm signal. When the ALARM SWITCH or BATTERY SWITCH is off, the alarm signal will not be activated.
  - \* N.B.: When you replace the used CO<sub>2</sub> cylinder, please use the CO<sub>2</sub> cylinder with Siphon Tube, not general CO<sub>2</sub> cylinder.
  - \* N.B: When the door is open during injection, the injection of CO<sub>2</sub> should be stopped automatically to protect the user

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### 6.3.3 How to set the temperature of the CO<sub>2</sub> backup system

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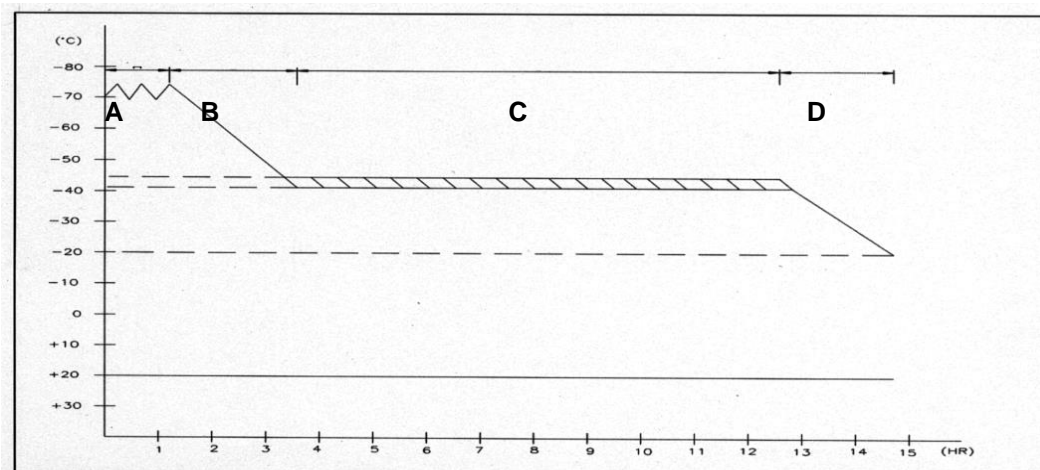
1. Push ① SET SWITCH
2. To change the set temperature, while you continue in pushing the button ① SET SWITCH, turn the ③ TEMP.ADJ. to the left and right to set the temperature,
3. Release the SET at the chosen temperature, the PRESENT lamp will be on and the present temperature of the chamber will be shown. The temperature is set to - 40 °C of factory value.

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### 6.3.4 Consuming time of the CO<sub>2</sub> gas (CO<sub>2</sub> gas 소모량)

**Tested by on the base of the following conditions: (다음 조건에 의한 테스트임)**

- Model: DF8517, 480 liter (모델 DF8517, 480 liter 기준)
- CO<sub>2</sub> cylinder capacity: 40 liter (CO<sub>2</sub> 봄베 용량: 40 liter 기준)
- CO<sub>2</sub> gas charge capacity: 40 liter (CO<sub>2</sub> Gas 충전량: 40 liter 기준)
- SET TEMP.: - 70 °C (설정 온도 : - 70 °C 기준)
- CO<sub>2</sub> backup system: SET TEMP. : - 40 °C (CO<sub>2</sub> backup system: 설정 온도: - 40 °C 기준)

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1. When the room temperature is above 30°C, the condenser cannot be cooled. It can cause malfunction of freezer.

=> If the room temperature around the unit is 28°C, please open the ventilation door and window to optimize the ventilation and cooling.

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2. When the room temperature is under 0°C, the oil of the compressor is condensed. It can have negative effect on the operation of the compressor. Each electrical part may have malfunction.

=> Under the circumstance of room temperature is below 0°C, make the place warm by using the other heating application and keep the temperature at least over 5°C.

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3. In case of the excessive dust around the unit, dust will pile up to the front of the condenser. This will cause the compressor to work inefficient, thereby causing compressor defect.

=> In case there is excessive dust around the unit, door which covers condenser should be cleaned up as often as monthly basis. Method of dust removal is wiping the dust off from the condenser by using soft brush.

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4. Under heavy humid circumstance, it can cause a technical malfunction for the electrical application and short circuit.

=> Ventilating the place properly by using the ventilator and opening the windows. An application, which can reduce moisture, will be more helpful.

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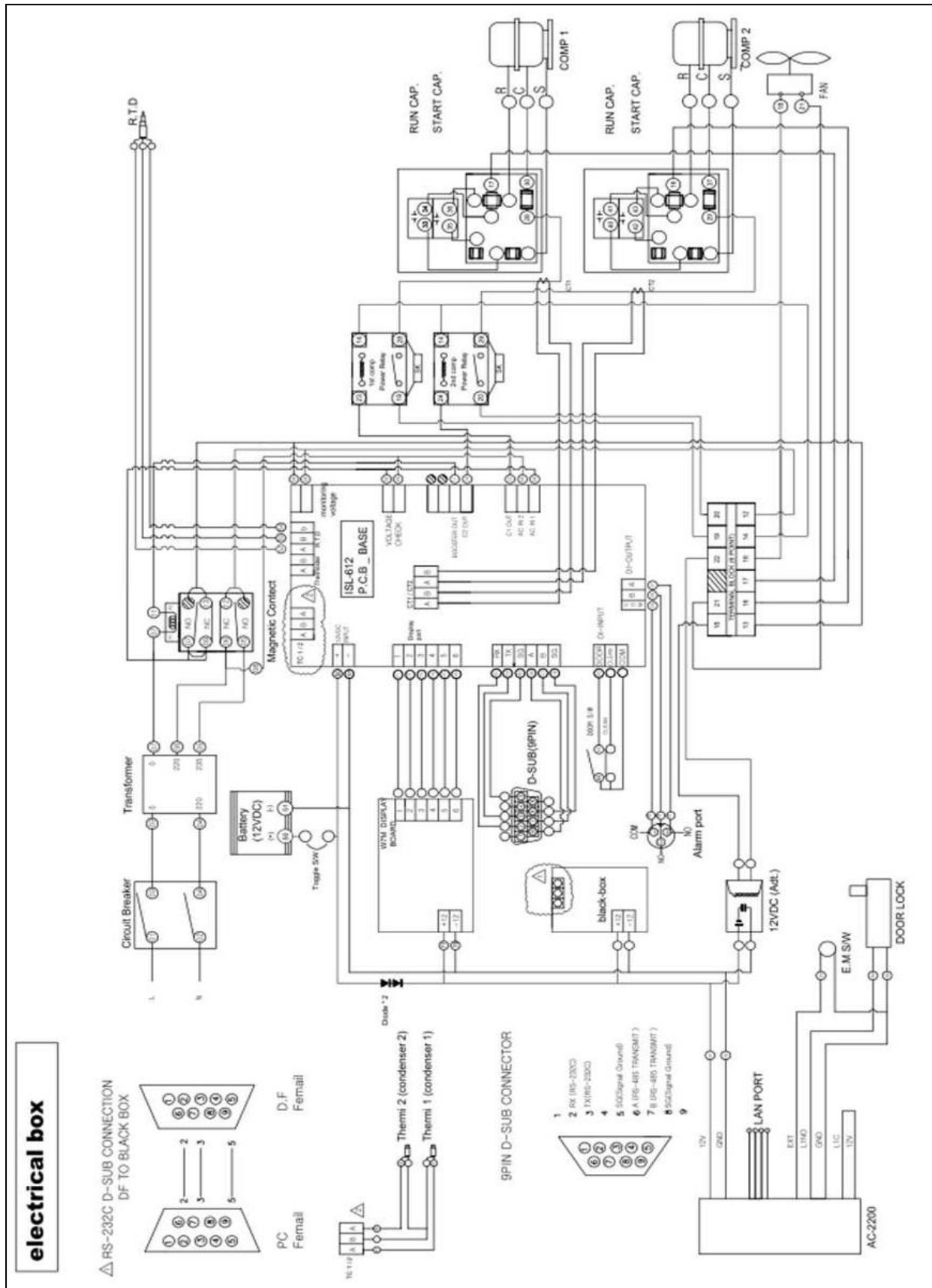
### 7.3 Electrical Schematic Diagram (전기 계통도)

## 7.4 Service Request (서비스 의뢰)

\* Before you register the service, please read the following information:

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### 1. The unit does not operate at all.

- Please check the power supply plug connection.
- Please check whether the temperature is set in a right way or not.
- Please check whether the unit is too close to a heating object or not.
- Please check whether the rear side of the unit is too close to the wall or not.
- Please check whether the condenser is blocked by dust or not  
(Remedy: Wipe off the dust by soft brush)
- Please check whether a big amount of test sample is inserted within a few hours or not.  
(Remedy: This is a temporal phenomenon, please wait)

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### 2. Running and abnormal noise

- Please check whether the floor is even or not.
- Please check whether the noise is coming from the compressor or from other source.
- \* The noise of the compressor is bigger than that of household refrigerators
- Do you hear a fan running sound?  
\* This unit is equipped with a fan motor which is used for air cooling system. When the fan motor is running, you can hear fan running noise.
- Do you hear a knocking sound(Creaking noise)?  
\* This sound originates from the parts of the unit, when they are expanded or shrunk.
- Do you hear water stream noise?  
\* The unit has refrigerant for it working fluid. You can hear circulation noise of refrigerant, which is not technical problem.

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## 3. A layer of ice has formed in the inside wall of the deep freezer

○ Put the samples with distances for good ventilation

○ Capping and sealing is required for samples with moisture.

○ Please checks door closing condition.

○ Please reduces the time of door open/close.

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## Warranty and contact information

### International Warranty

The warranty period starts on the moment the unit leaves the production facility. After this date one year warranty on the parts is valid. During this first year component parts proven to be defective could be send to your local dealer and will be covered by the first year warranty on parts only. Expendable items, i.e., glass, filters, pilot lights, light bulbs and door gaskets are excluded from this warranty. ilShinbiobase Co. Ltd. shall not be liable for any indirect or consequential damages including, without limitation, damages relating to lost profits or loss of products.

### Conditions and terms of sale

All prices are ex-works from the ilShinbiobase Co. Ltd. Factory in Korea, excluding insurance, seaworthy packing and customs duty. Effective prices are those valid on date of confirmation of order. All specifications mentioned in the price list are based on 230 V, 50/60Hz power supply. Continuing research and improvements may result in specification changes. Errors and misprints accepted.

### Contact information

In case service of sales back up is needed please do not hesitate to contact your local ilShin Dealer:

--

Or contact the ilShinbiobase Co. Ltd. Headquarter:

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