# Pocket Freezer Series

Users Manual, Ver. 2.0

04/04/17

**Failure to read, understand and follow the instructions in this manual may result in damage to the unit, injury to operating personnel and poor equipment performance.**

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These operating instructions are protected by copyright. Rights resulting thereof, particularly reprint, Photomechanical or digital post processing or reproduction, even in part, are only allowed with the written consent of ilShinbiobase.

This regulation does not apply to reproductions for in-plant use.

### **Trademarks**

All trademarks mentioned in the operating instructions are the exclusive property of the respective

ilShinbiobase Co,Ltd. provides this document to its customers with a product purchase to use in the product operation. This document is copyright protected and any reproduction of the whole or any part of this document is strictly prohibited, except with the written authorization of ilShinbiobase Co,Ltd.

The contents of this document are subject to change without notice. All technical information in this document is for reference purposes only. System configurations and specifications in this document supersede all previous information received by the purchaser.

ilShinbiobase Co,Ltd. makes no representations that this document is complete, accurate or error-free and assumes no responsibility and will not be liable for any errors, omissions, damage or loss that might result from any use of this document, even if the information in the document is followed properly.

This document is not part of any sales contract between ilShinbiobase Co,Ltd. and a purchaser. This document shall in no way govern or modify any Terms and Conditions of Sale, which Terms and Conditions of Sale shall govern all conflicting information between the two documents.

**© ilShinbiobase Co,Ltd. All rights reserved.**

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**(Read and understand the manual before using this machine)**

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**(While machine works, the printed sign can be detached by bad sticking then, attach the same one)**

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**Only authorized personnel may service this equipment**

**Before Installation**

## -Transportation-

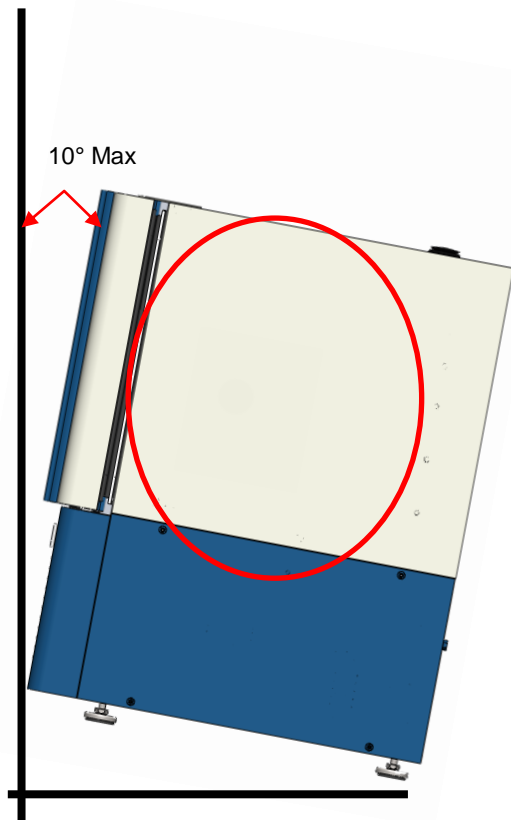
If the unit is tilted excessively, the compressor lubricant may flow from the compressor to the line system. An insufficient filling level may damage the compressor when the refrigeration system is started.

**Transport the unit only vertical position.** When transporting, the unit must not tilt beyond 10° determined.

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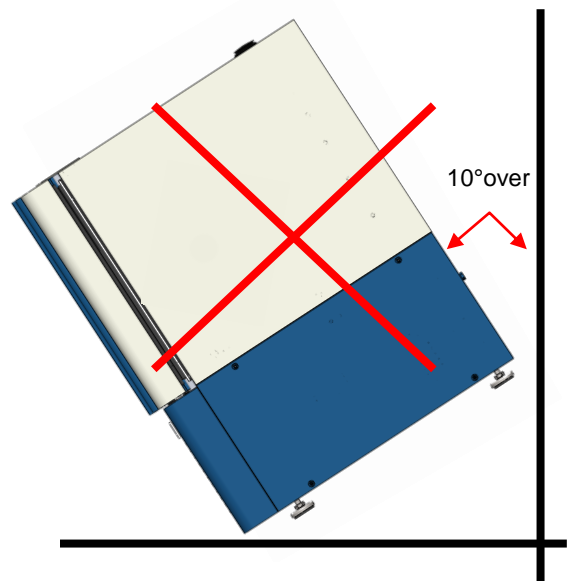
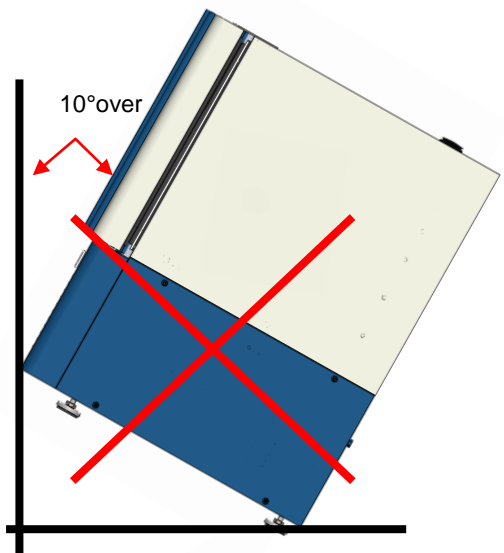
- Recommended room temperature: between 5℃ and 28 ℃. If the room temperature is normally > 30 ℃ please ensure that suitable air-conditioning is installed.
- Make sure the floor is level and even.
- Make sure the condenser grid is not blocked. (Adequate air circulation is required)
- Install the two spacer pins on the back of the freezer to allow heat dispersion behind the freezer. (fig.1)

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### **2. Turning Freezer on(Freezer 켜기)**

- Push the power plug firmly into the power socket. (Please check you have the correct voltage)
- Make sure the main power switch at the back of the freezer is switched on. (Switched upwards) (fig.2)

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### **3. Recommended set points(권장 설정온도)**

- a. set-point – 80 °C
- b. These set points are factory set and gives the freezer the optimum conditions for prolonged life

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**We wish you a quick and trouble free installation!**

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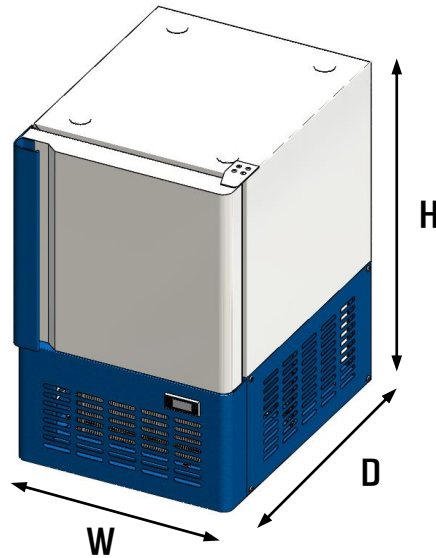
---

## 1. Dimension of The Freezer Models (모델 별 규격)

### 1.1 layout

Model No.	Exterior Dimensions	Interior Dimensions	Capacity(L)
	WxDxH	WxDxH	
DF8502	470x620x780mm	300x440x300mm	40
DF8503	760x620x780mm	600x440x300mm	80

The operating temperature of Pocket Freezer is  $-65^{\circ}\text{C}$  to  $-86^{\circ}\text{C}$  (normal temperature for usages is  $-70^{\circ}\text{C}$  to  $-80^{\circ}\text{C}$ ). In addition, if the last two numbers(<sup>xx</sup>) of the model are the same, the same operating temperature is applicable.

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## 2. Specification(기술 사양)

### 2.1 Deep Freezers Characteristic (초저온 냉동고 특징)

1. Hermetically sealed cooling system with very durable compressors
2. Efficient evaporator with maximum contact with walls for fast cooling down
3. polyurethane foamed insulation with VIP© insulation
4. Fixed support with built in adjustable stands ( unique )
5. Cabinet with round corners for easy cleaning
6. 2 Built-in access ports
7. Very user-friendly control panel
8. Easy to clean condenser filter
9. Quiet running fans and compressors create a very low noise level (50~65 dbA)

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### 2.2 Specification

Model No.	Capacity (L)	Weight (Kg)	Low Temp. °C	Usage Temp. °C	Electrical (V, Hz)	Power (Hp)	Insulation	Const- ruction
DF8502	40	79	-86	-60 ~ -80	220/230, 50/60	3/8	High Density Polyuretha ne With VIP	Textured steel or stainless steel (optional)
DF8503	80	99				5/8		

### 3. Installation (설치)

#### 3.1 Installation place (설치 장소)

1. Not to subject to direct sunlight
2. Adequate air circulation
3. Not subject to excessive dust
4. Not subject to mechanical vibration or shock
5. Ambient temperature: 5℃ to 33℃, Optimum temp. is 18℃ to 25℃
6. Operating Height : below 2,000m
7. Operating moisture: below 80% RH, in case that the max. operating temperature is 33℃, it should be below 57% RH
8. Allowance of supplying voltage change : lower than 10%

\* Caution: In case that you install the unit other places than mentioned, the unit cannot be operated properly.

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#### 3.2 How to install (설치 방법)

1. Keep the distance 15cm or longer between the wall and the backside of the unit.
2. Do not install the unit in an uneven position.
3. The ventilation should not be obstructed

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## 4. Power supply, how to earth and main power switch

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1. To ensure trouble-free connection, connect the plug firmly in the power socket.
2. Don't connect the plug of the unit together with plug of other units in the same socket.
3. Make sure that the power supplying cord is not folded and pressed.
4. If you use an extension wire, it should has capacity above 250V, 15A and not longer than 3m.
5. This unit is exclusively designed for 220(230)V, 60(50)Hz, please check the operating voltage, before use. Under the 110V, transformer for elevating voltage (above 3Kw) should be used.

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#### 4.2 How to earth (접지 방법)

1. For security reasons, the unit should be earthed.
2. If the power socket is fitted with earth rod (230V), please check the outlet for earth before use.
3. If the power socket is not fitted with earth rod, first, connect the earth wire to the unit. Then, connect it to a bronze plate. Subsequently, bury it in the moistened ground. (25 cm under the ground or deeper)

\* Warning: Do not earth to a gas pipe, a telephone line or lightning conductor.

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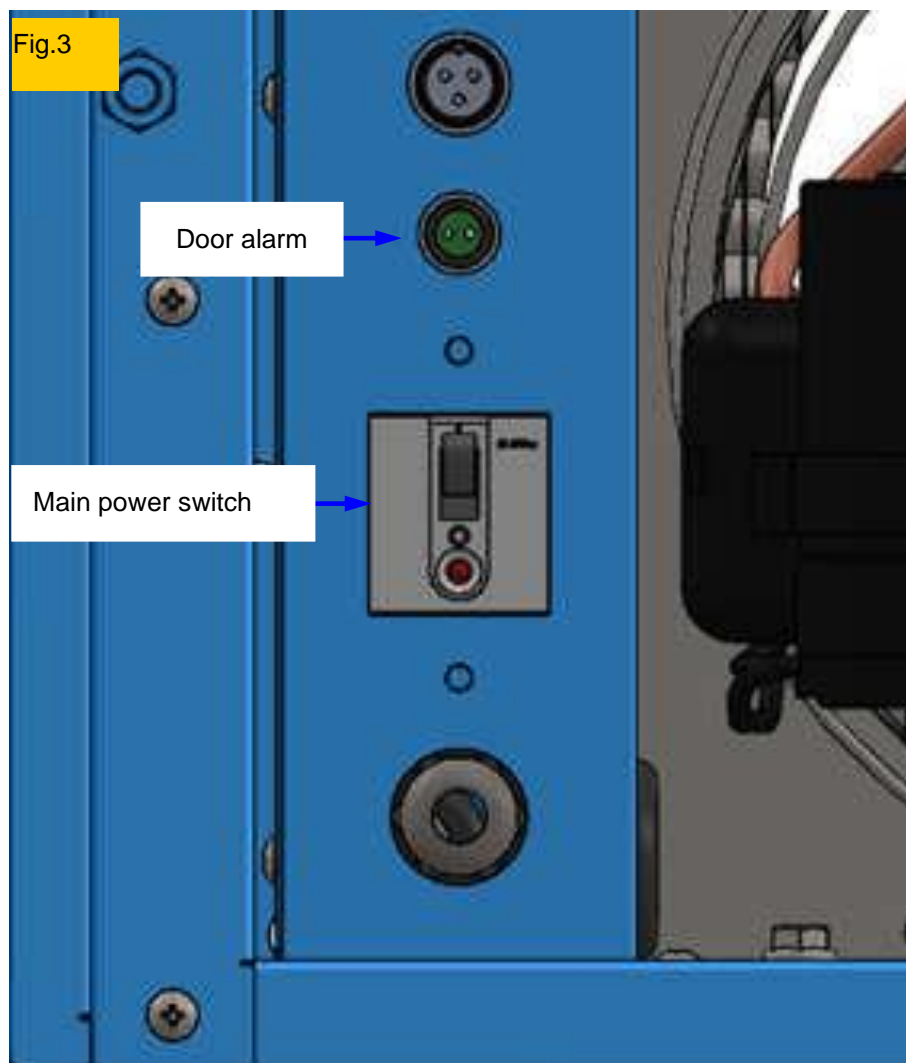
The main power switch is located rear the product next to the compressor. After power supply is checked and the unit is earthen. By switching the main power switch in **ON** position the Freezer will use the power and on this time is ready for operation. If the freezer is not in use, please put the main switch in **OFF** position.

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#### 4.4 Door Alarm(도어 알람)

The door alarm can be used in combination with the CO<sub>2</sub> backup system. It is a safety to make sure the operator will not get liquid CO<sub>2</sub>/LN<sub>2</sub> in the face in case of opening the door of the freezer. Always connect the CO<sub>2</sub>/LN<sub>2</sub> backup system to the door alarm connection on the back of the freezer.

For safety measure resistance on the contacts with Ohm meter

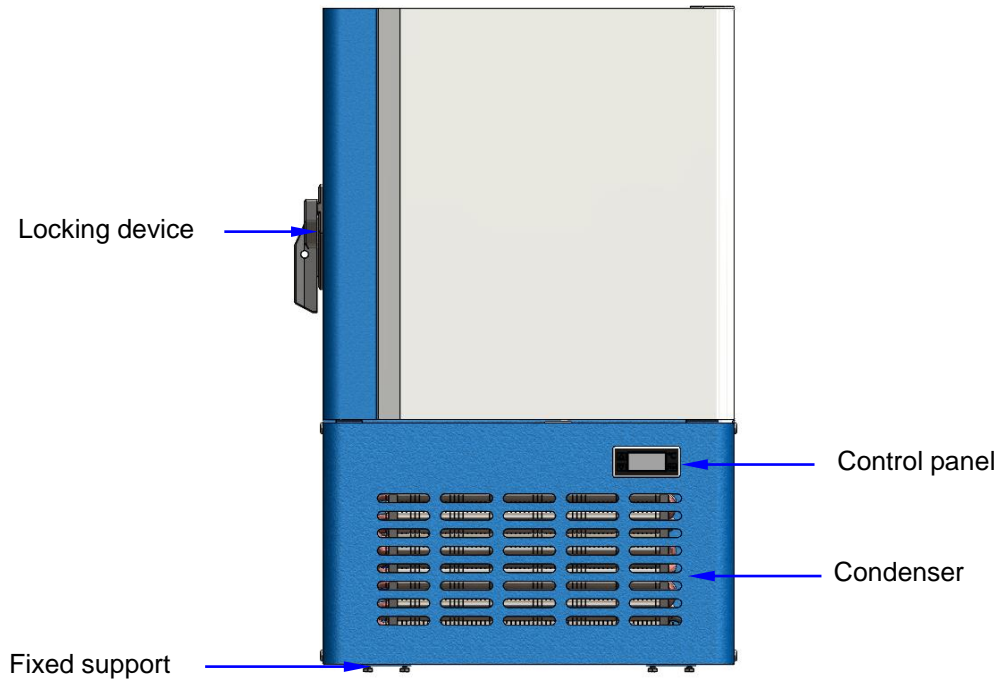
1. do contact 1&2 show resistance? < 0 Ohm> It means Door is Closed.
2. open the door (chamber)
3. do contact 1&2 show unlimited resistance? It means Door is Opened.

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

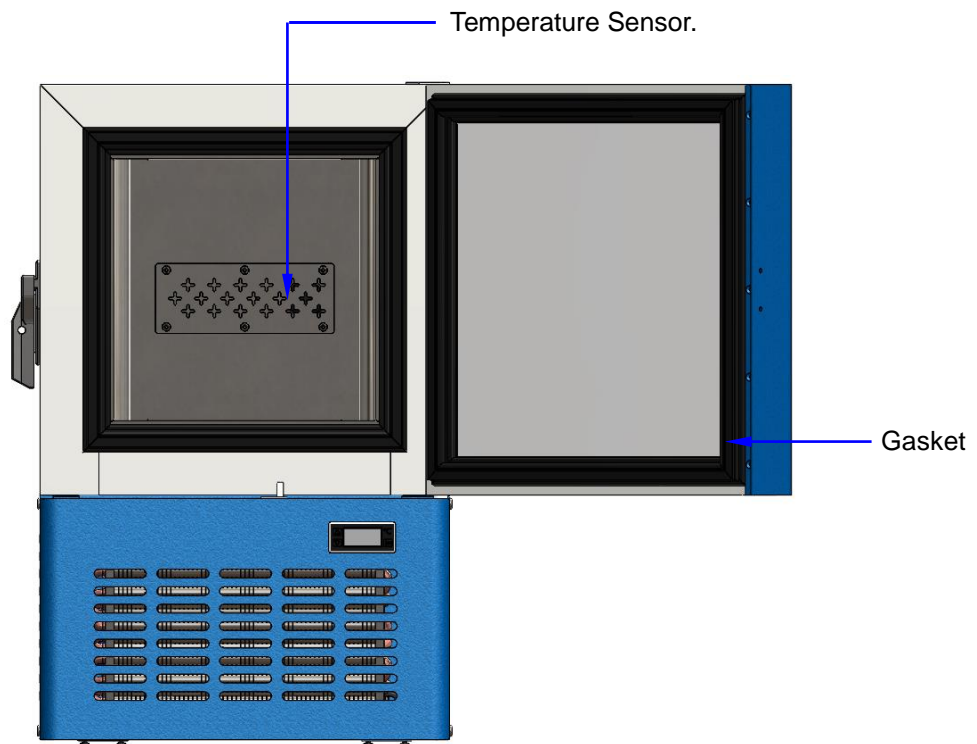
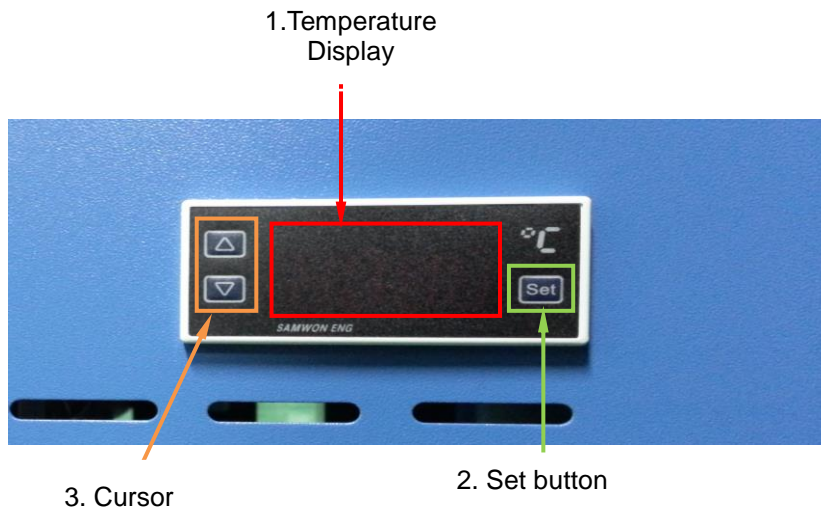
## 5.6 How to Operate (작동 순서)

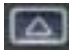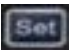1. Turn on the circuit breaker. (See the rear guard)
2. All functions will be operational. The signal will be on and the inside temperature of the Chamber will be shown.
3. Push the set button  once to check the temperature. The SET POINT will be shown. (To adjust SET POINT, use the CURSOR button  )
4. Press the SET button  again to end the setting.

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### 6.2 Fireboard Storage Boxes and Dividers(박스&칸막이)

#### **External dimensions:**

External dimensions of 2 inch boxes (model PP0060): 133x133x49 mm (DxWxH)

External dimensions of 3 inch boxes (model PP0061): 133x133x77 mm (DxWxH)

#### **Internal Dimensions:**

Internal dimensions of 2 inch boxes (model PP0060): 124x124x45 mm (DxWxH)

Internal dimensions of 3 inch boxes (model PP0060): 124x124x70 mm (DxWxH)

## Possible dividers:

Dividers are also made of plasticized card box. Height of all dividers is 35 mm

### **Model: Description:**

PP1049 Storage box divider 7 x 7

PP1064 Storage box divider 8 x 8

PP1081 Storage box divider 9 x 9

PP1100 Storage box divider 10 x 10

Thick Plasticized Card box Material

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1. When the room temperature is above + 30°C. When the room Temperature is above +30°C, the condenser cannot be cooled. It can be one of cause of trouble.

=> If the room temperature around the unit is +28°C, please open the condenser door and window to optimize the ventilation and cooling.

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2. When the room temperature is under 0°C. The oil of the compressor is condensed. It can have negative effect on the operation of the compressor. Each electrical part can be wrongly operated

=> Under the circumstance that the room temperature is under 0°C, make the place warm by using the other heating application and keep the temperature minimal +5°C.

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3. In case there is excessive dust around the unit. This unit use the condenser which is air cooled, In case of the excessive dust around the unit, dust will pill up the front of the condenser. This will cause the compressor to nor work efficiently, This is one of the cause of a defect compressor

=> In case there is excessive dust around the unit door which covers condenser should be often once in a month and remove dust and clean up. Also wipe the dust off the condenser with a soft brush.

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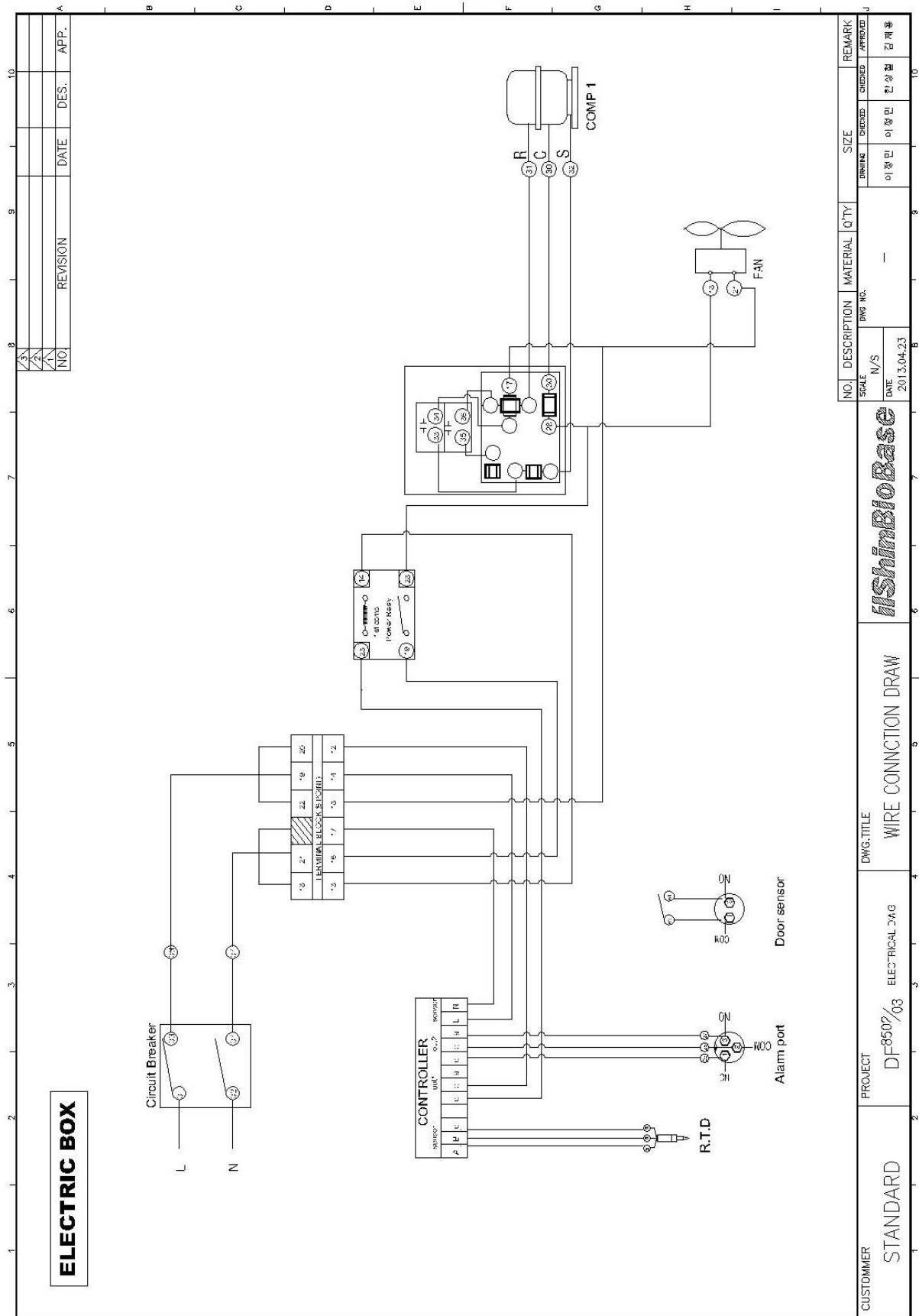
4. Under heavy humid circumstances. It can cause a technical malfunction in the electrical application and short circuit.

=> Ventilation the place properly by using the ventilator and opening the windows. An application, which can reduce moisture, will be helpful.

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## 7.2 Electrical Schematic

## 7.3 Registration of Service (서비스 의뢰)

\* Before you register for the service, please read the following information:

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### 1. The unit does not operate at all.

○ Please, checks the power supply (230V) plug connection.

○ Please, checks whether the temperature is set in a right way.

○ Please, checks whether the unit is too close to a heating object.

○ Please, checks whether the rear side of the unit is too close to the wall.

○ Please, check whether the condenser is blocked by dust

(Remedy: wipe off the dust by soft brush)

○ Please, checks whether a big amount of test sample is inserted within a few hours. (Remedy: this is a temporal phenomenon, please just wait)

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### 2. Loud running noise

○ Please, check whether the floor is even.

○ Please, check whether the noise is coming from the compressor.

\* The noise of the compressor is bigger than that of household refrigerators

○ Do you hear a turbo sound?

\* This unit is equipped with a fan motor which is used for air cooling system. When the fan motor is running, you can hear some noise.

○ Do you hear a knocking sound?

\* This sound originates from the parts of the unit, when they are expanded or shrunk.

○ Do you hear water stream?

\* The unit is equipped with circulation of refrigerant. You can hear the refrigerant stream through the circulation. It is not a technical problem.

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## 3. A layer of ice has formed in the inside wall of the deep freezer

○ Put the samples with distances for good ventilation

○ If a sample contains much moisture, then put a top on it, or keep it in wrap.

○ Please, checks whether the door is closed.

○ Please reduces the times to open the door.

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### Warranty and contact information

#### International Warranty

The warranty period starts on the moment the unit leaves the production facility. After this date one year warranty on the parts is valid. During this first year component parts proven to be defective could be send to your local dealer and will be covered by the first year warranty on parts only. Expendable items, i.e., glass, filters, pilot lights, light bulbs and door gaskets are excluded from this warranty. IlShinbiobase Co., Ltd shall not be liable for any indirect or consequential damages including, without limitation, damages relating to lost profits or loss of products.

#### Conditions and terms of sale

All prices are ex works from the ilShinbiobase Co. Ltd Factory in Korea, excluding insurance, seaworthy packing and customs duty. Effective prices are those valid on date of confirmation of order. All specifications mentioned in the price list are based on 230 V, 50/60Hz power supply. Continuing research and improvements may result in specification changes. Errors and misprints accepted

#### Contact information

In case service of sales back up is needed please do not hesitate to contact your local ilShin Dealer:

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Or contact the ilShinbiobase Co., Ltd Headquarter: