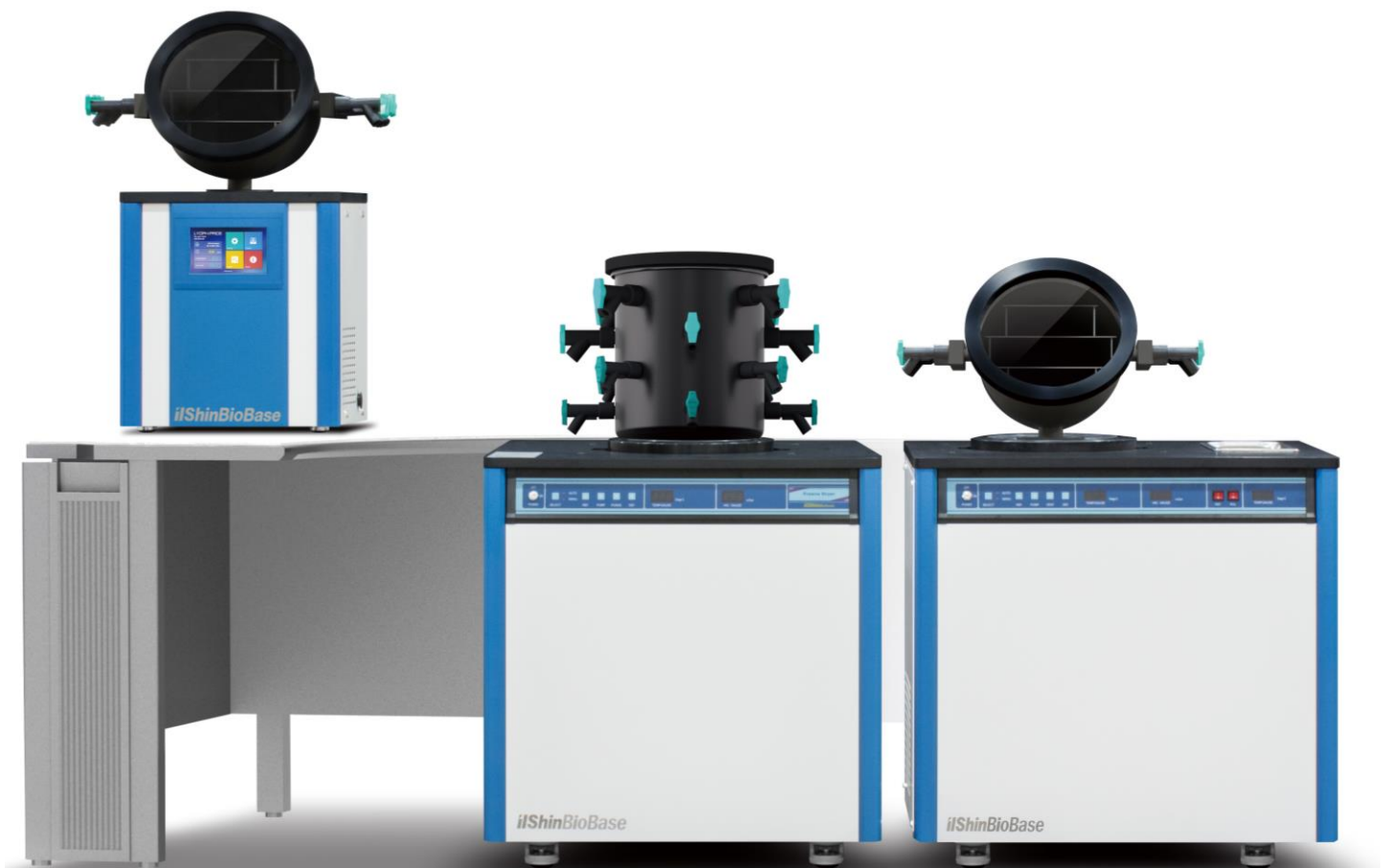
Freeze Dryer Series

Users Manual, Ver 4.0

03/04/17

Failure to read, understand and follow the instructions in this manual may result in damage to the unit, injury to operating personnel and poor equipment performance.

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These operating instructions are protected by copyright. Rights resulting thereof, particularly reprint, photomechanical or digital post processing or reproduction, even in part, are only allowed with the written consent of ilShinbiobase.

This regulation does not apply to reproductions for in-plant use.

Trademarks

All trademarks mentioned in the operating instructions are the exclusive property of the respective.

ilShinbiobase Co. Ltd. provides this document to its customers with a product purchase to use in the product operation. This document is copyright protected and any reproduction of the whole or any part of this document is strictly prohibited, except with the written authorization of ilShinbiobase Co. Ltd.

The contents of this document are subject to change without notice. All technical information in this document is for reference purposes only. System configurations and specifications in this document supersede all previous information received by the purchaser.

ilShinbiobase Co. Ltd. makes no representations that this document is complete, accurate or error-free and assumes no responsibility and will not be liable for any errors, omissions, damage or loss that might result from any use of this document, even if the information in the document is followed properly.

This document is not part of any sales contract between ilShinbiobase Co. Ltd. and a purchaser. This document shall in no way govern or modify any Terms and Conditions of Sale, which Terms and Conditions of Sale shall govern all conflicting information between the two documents.

© ilShinbiobase Co. Ltd. All rights reserved.

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(Read and understand the manual before using this machine)

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(While machine works, the printed sign can be detached with poor adhesion, then attach with the same one)

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Only authorized personnel may service this equipment

Before Installation

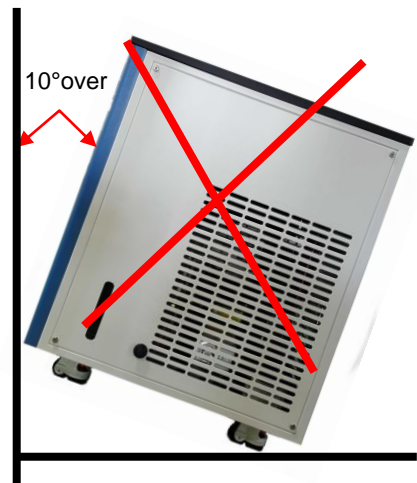
-Transportation-

If the unit is tilted excessively, the compressor lubricant may flow from the compressor (or oil separator) to the line system. Insufficient filling level may damage the compressor when the refrigeration system is started. Keep vertical position while transporting the unit. During transportation, the unit must not tilt more than 10° .

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
Max. tilt angle transportation (Enable position)

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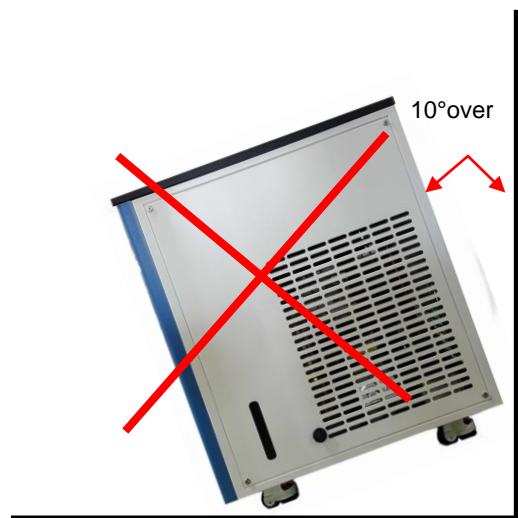1. Recommended room condition (권장 장소)

- a. Recommended room temperature: between 5℃ and 28℃. If the room temperature is over 30℃, please ensure that suitable air-conditioning is provided.
- b. Make sure the floor is level and even.
- c. Make sure the condenser grille is not blocked. (Adequate air circulation is required)

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2. Turning Freeze Dryer on(Freeze Dryer 켜기)

- a. Push the power plug firmly into the power socket.(Please check you have the correct voltage.)
- b. Turn the Key switch to the right (clockwise) 90 degrees and set it to ON.
- c. Please check your monitor

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- a. We recommend the condenser grille is free from dust every month.
- b. We recommend a maintenance visit by your local dealer every year.

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We wish you a quick and trouble-free installation!

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1. Specifications of the model(모델별 사양)

A. FD8508 – FD8518

		FD8508	FD8512	FD8518
Ice capacity		5 L	10 L	12 L
Totally Ice capacity		8 L	12 L	18 L
Condenser Temperature		-70 to -85 °C		
Ext. Dimensions		800 x 705 x 860 mm		
Refrigeration		1 1/4 HP		
Electrical		220V 60Hz, 230V 50Hz		
Option	Drying Chamber	Multi Drying Chamber / Drying Chamber / Clear Drying Chamber / Heated Drying Chamber / Manifold Chamber		
	Product Shelf	PH1316		
	Vacuum pump	ISL136(196LPM) / ISL201(283LPM)		

B. FDS8508 – FDS8518

		FD8508	FD8512	FD8518
Ice capacity		5 L	10 L	12 L
Totally Ice capacity		8 L	12 L	18 L
Condenser Temperature		-70 to -85 °C / Shell Freezer : -40 °C		
Ext. Dimensions		850 x 750 x 860 mm		
Refrigeration		1 1/4 HP + 3/8 HP		
Electrical		220V 60Hz, 230V 50Hz		
Option	Drying Chamber	Multi Drying Chamber / Drying Chamber / Clear Drying Chamber / Heated Drying Chamber / Manifold Chamber		
	Product Shelf	PH1316		
	Vacuum pump	ISL136(196LPM) / ISL201(283LPM)		

**C. MCFD8508
– MCFD8518**

		FD8508	FD8512	FD8518
Ice capacity		5 L	10 L	12 L
Totally Ice capacity		8 L	12 L	18 L
Condenser Temperature		-70 to -85 °C		
Ext. Dimensions		1005 x 750 x 885 mm		
Refrigeration		1 1/4 HP		
Electrical		220V 60Hz, 230V 50Hz		
Option	Drying Chamber	Multi Drying Chamber / Drying Chamber / Clear Drying Chamber / Heated Drying Chamber / Manifold Chamber		
	Product Shelf	PH1316		
	Vacuum pump	ISL136(196LPM) / ISL201(283LPM)		

**D. TFD8501
- TFD8503**

		TFD8501	TFD8503
Ice capacity		1 L	2 L
Totally Ice capacity		2 L	3 L
Condenser Temperature		-70 to -85 °C	
Ext. Dimensions		450 x 525 x 500 mm	800 x 650 x 400 mm
Refrigeration		3/8 HP	
Electrical		220V 60Hz, 230V 50Hz	
Option	Drying Chamber	Multi Drying Chamber / Drying Chamber / Clear Drying Chamber / Heated Drying Chamber / Manifold Chamber	
	Product Shelf	PH1212	
	Vacuum pump	ISL6(100LPM) / ISL136(196LPM)	

2. Installaion (설치)

2.1 Installation place (설치장소)

1. Not to subject to direct sunlight
2. Adequate air circulation
3. Not subject to excessive dust
4. Not subject to mechanical vibration or shock
5. Operating temperature: + 5 °C to + 33 °C
6. Operating height: below 2.000 m
7. Operating moisture: below 80% (in case that the max. operating temperature is + 33 °C , it should be below 57 %)
8. Allowance of supplying voltage change: lower than +/- 10%

*** In case that you install the unit at any other places than mentioned, the unit Can't operate properly.**

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2.2 How to install (설치방법)

1. Keep a distance of 30 cm or longer between the wall and the backside of the unit.
2. Do not install the unit at an uneven position.
3. Make the unit fixed with fixing screws on a steep floor.
4. The ventilation should not be obstructed.

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1. To ensure trouble-free connection, connect the plug firmly in the power socket.
2. Don't connect the plug of the unit together with plug of other units in the same socket.
3. Make sure that the power supplying cord is not folded and pressed.
4. If you use an extension wire, it should have a capacity above 230V, 15A and not longer than 3 m.
5. This unit is exclusively designed for 230V, 50Hz, please check the operating voltage, before use.

*** Under the 110V, the transmitter for elevating voltage (above 3Kw) should be used**

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4.1 Index of the parts for the FD8508 ~ FD8518 (부분별 명칭)

4.2 Index of the parts for the FDS8508 ~ FDS8518 (부분별 명칭)

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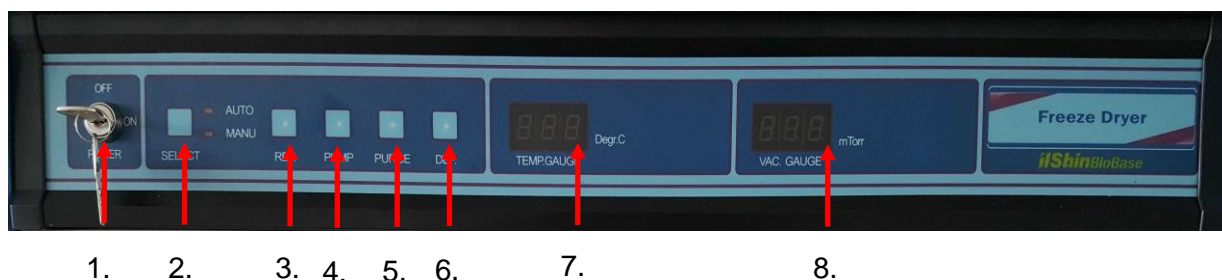
5.2. Manual Mode (수동 모드)

Before drying of the test sample

1. Turn the Key switch to the right (clockwise) 90 degrees and set it to ON.
2. Now, you can operate all functions. The MANU lamp will be on and the present Temperature of the cold trap will be indicated on the TEMP. GAUGE.

*** No value will be indicated on VAC. GAUGE on this time.**

3. Press (3) REF. Button once, the lamp will be on and the refrigeration of the unit will be operating.
4. After about 10 min, the number of the TEMP GAUGE will go down to freezing point.
5. If the temperature is under the -40°C , press the (4) PUMP button once.
6. The lamp will be on and the vacuum pump will be operating.
7. After about 10 minutes, the value "999" will show on the Vac. Gauge and this value will go down to approximately 10mTorr .
8. From this point, the test sample will be dried gradually.

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When the test sample is dried:

1. When the test sample is dried, push the button (3) REF. Once. The REF led will go out and the refrigeration will stop.
 2. Push the button (4) PUMP button, then the PUMP led will go out and the vacuum pump will stop.
 3. Push the (5) VENT button once, then the VENT led will go out and the vacuum state of the chamber and vacuum valves will be released by the automatic valve. (During the switch-off of the vacuum, little noisy sound can be heard)
 4. After the system is totally released, push the button (5) VENT once more to go to the next drying process. The VENT led will go out and the automatic valve will be closed.
 5. Push the (6) DEF. Button once, then, the DEF led will go on and the ice on the cold trap will be defrosted by the Hot Gas system.
 6. After the ice is totally defrosted, push the (6) DEF button once more. The DEF led will go out and the heating process around the chamber will stop.
- * If the next drying process is not directly followed, you could better defrost the ice in the cold trap without usage of the DEF button

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5.3 Automatic Mode (자동 모드)

Before drying of the test sample:

1. Turn the key switch to the right (clockwise) 90 degrees and set it to ON.
2. Now, you can operate all functions. The MANU lamp will be on and the present temperature of cold trap will be indicated on the TEMP. GAUGE.
* No value will be indicated on VAC. GAUGE on this time
3. Push the (2) SELECT button once, then the MANU led will be turned off and the AUTO led will be on. The freezer will be operated, while the led of the (3) REF. button is on.
4. After about 10 min, the value of the TEMP GAUGE will go down to freezing point. If the temperature is under the -40°C , the (4) PUMP led will be on by the automatic sensor and the vacuum pump will be operated
5. After about another 10 minutes, a value of "999" will be shown in the VAC. GAUGE display and this value will slowly go down to approximately 10mTorr
6. From this point, the test sample will be dried gradually.

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When the test sample is dried:

1. When that the test material is dried, push the button (2) SELECT once. The AUTO led will be turned off and The MANU led will be on.
The led of (3) REF. Button and the (4) PUMP button will be turned off and all functions will stop.
 2. Push the button (4) PUMP button, then the PUMP led will go out and the vacuum pump will stop.
 3. Push the (5) VENT button once, then the VENT led will go on and the vacuum state of the chamber and vacuum valves will be released by the automatic valve. (During the switch-off of the vacuum, little noisy sound can be heard)
 4. After the system is totally released, push the button (5) VENT once more to go to the next drying process. The VENT led will go out and the automatic valve will be closed.
 5. Push the (6) DEF. Button once, then, the DEF led will go on and the ice on the cold trap will be defrosted by the Hot Gas system.
 6. After the ice is totally defrosted, push the (6) DEF button once more. The DEF led will go out and the heating process around the chamber will stop.
- * If the next drying process is not directly followed, you could better defrost the ice in the cold trap without usage of the DEF button

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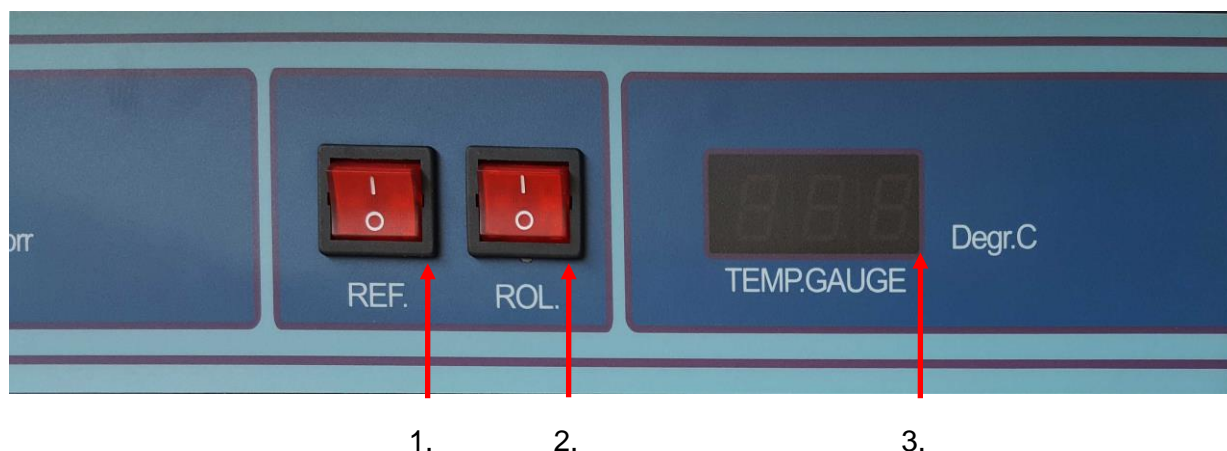
5.4 Shell freezer operation (Shell Freezer 작동법)

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
5.4.2. How to operate the shell freezer (작동 방법)

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5.5 Freeze Dryer With Micro Concentrator operation (MCFD작동법)

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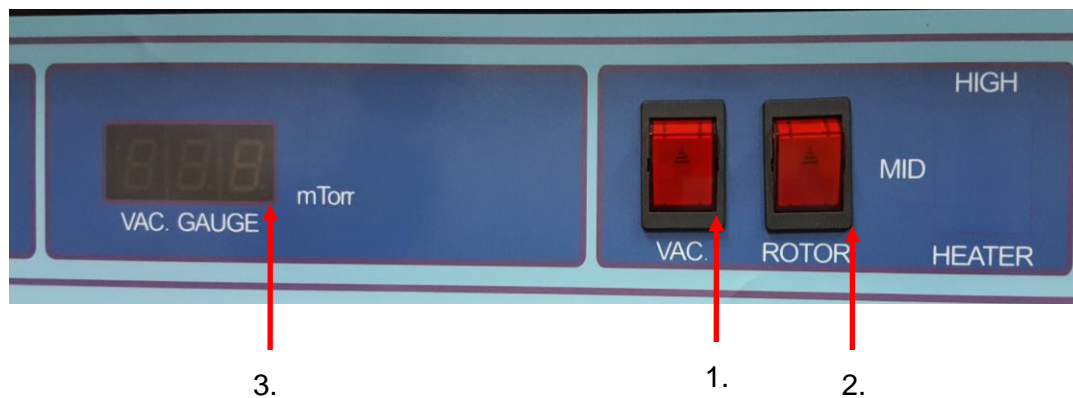5.5.2. How to operate the shell freezer (작동 방법)

- Now, you can operate all functions. Push the vacuum button(1) once activates the Vacuum of the Micro Concentrator, Vacuum pressure appears on the gauge(3).
 - Push the ROTOR button(2) is pressed once, the Micro Concentrator operates (rotates).
- 진공 버튼(1)을 누르면 Micro Concentrator의 Vacuum이 활성화되고, 진공 압력이 게이지 (4)에 나타납니다.
 - ROTOR 버튼을 1회 누르면, Micro Concentrator가 동작(회전)합니다.

6. Accessories(악세서리)

6.1 Configuration (구성)

Accessories Specifications.

1) Multi Drying Chamber

Model	Discription
MDC 1208	8 x ½" Valves, 275(H) x 230mm(D), 304 Stainless steel, 4 Shelves
MDC 1208 H	MDC 1208 with heating shelves (+40°C at ambient Temp.)
MDC 1208 HS	MDC 1208 H with stoppering system

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3) Clear Drying Chamber


Model	Discription
CDC 3332	2heated shelves 330mm(H) x 330mm Dia
CDC 3333	Stoppering and 2heated shelves 330mm(H) x 330mm Dia

4) Heated Drying Chamber

Model	Discription
HDC 1212	12 x ½" Valves, 275(H) x 230mm(D), 304 Stainless steel
HDC 1316	8 x ½" and 8 x ¾" Valves 320 mm (H) x 300 mm (D) with Heater function
HDC 1218	18 x ½" Valves 320 mm (H) x 300 mm (D) with Heater function

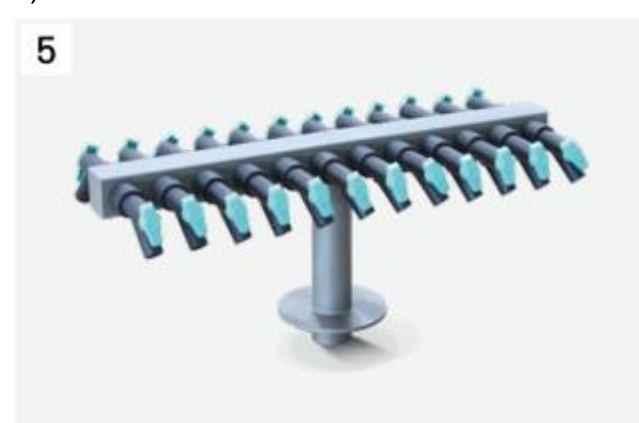5) Manifold Chamber

Model	Discription
MC 1224	12 x ½" Valves 275mm (H) x 230 mm (D) with Heater function
MC 3412	12 x ¾" Valves
MC 1324	12 x ½" and 12 x ¾" Valves
MC 1204	4 x ½" Valves

6) Product Heater

Model	Discription
PS1212	3 Shelves for DC 1212
PS1316	3 Shelves for DC 1316/1218
PHS1212	3 Heated Shelves for DC 1212
PH1316	3 Heated Shelves for DC 1316/1218


7) Specifications of Flask

Size	Complete	Bottom	Top	Diameter	Bottom H/I.D.
40ml	CF0040	FB0040	FT0034	1/2"	76 x 34(mm)
80ml	CF0080	FB0080			115 x 34(mm)
120ml	CF0120	FB0120	FT0059	3/4"	69 x 60(mm)
150ml	CF0150	FB0150			85 x 60(mm)
300ml	CF0300	FB0300			145 x 60(mm)
600ml	CF0600	FB0600	FT0090		135 x 90(mm)
900ml	CF0900	FB0900			190 x 90(mm)
1200ml	CF1200	FB1200			240 x 90(mm)

6.2. Vacuum pump(진공펌프)

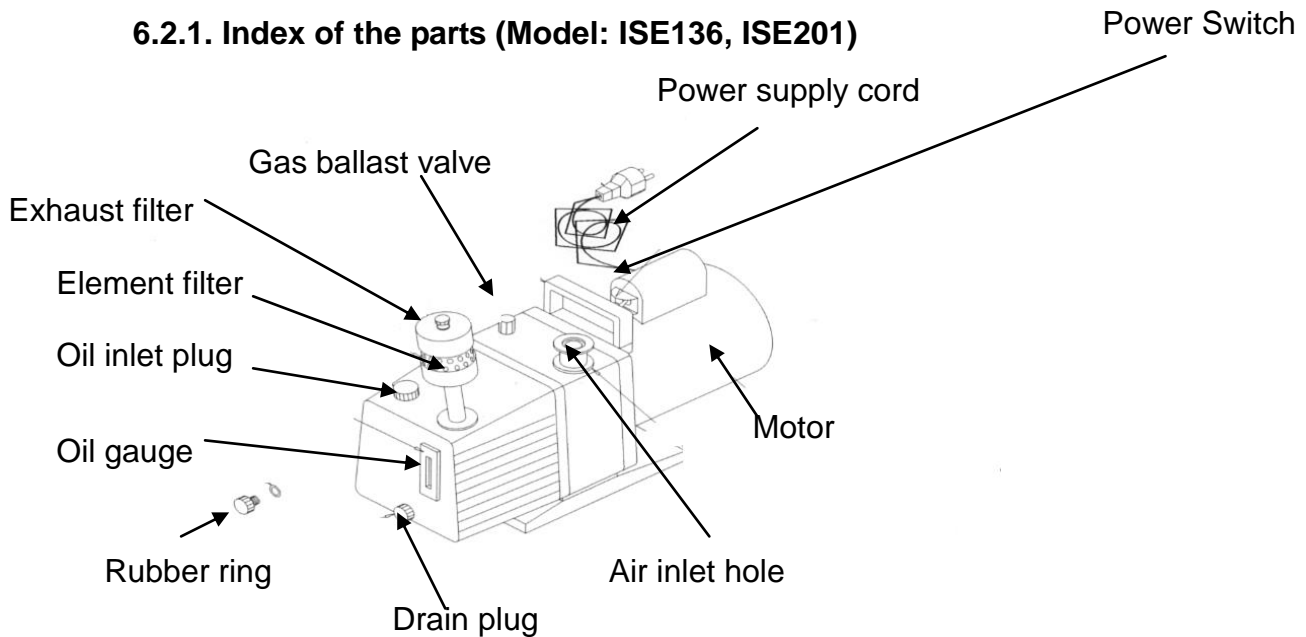6.2.1. Index of the parts (Model: ISE136, ISE201)

6.2.2. Oil change frequency and how to change the oil(오일교환시기 및 방법)

- The oil of the vacuum pump should be changed over a certain period to make sure the pump will run optimally (just like the engine oil of a car should be changed after a certain km)
- The oil should be changed when the color of the oil seems to be not clear anymore or dark yellow to brown tint through the oil gauge.
- The changing period of the oil depends on the test sample which is freeze dried, but a general rule of thumb is to change the oil after every 700 to 1.000 running hours.

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How to change the vacuum pump oil (오일 교환 방법)

Step 1: After making the clasper connected to the vacuum tube loose, disconnect the pump from the vacuum circuit.

Step 2: Put a basin under the oil drain plug. Make the drain plug open while turning the plug to the left (counter clockwise). Please, check whether the oil is fully drained by checking the oil gauge.

* Do not lose the rubber ring in the front part of the drain plug.

Step 3: Please, check the rubber ring in the front part of the plug. If it is not present, please, insert it precisely.

Step 4: Turn the oil drain plug to the right (clockwise) firmly.

* N.B. the drain plug and pump drain should be fit on the right way.

Step 5: Turn the filling plug of the pump to the left (counter clockwise) to open. Fill the new oil up through the opening.

Step 6: Fill up to the maximum line of oil gauge (For Model VP8915, VP8920: 1.3 liter)

Step 7: If the filling up is finished, please turn the filling plug to the right (clockwise) and close it firmly.

* N.B. the drain plug and pump drain should be fit on the right way.

Step 8: Bring the vacuum pump back and connect the vacuum line of the unit to the pump precisely with the clamp.

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6.2.3 How to change elements of exhaust filter

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- When the solvent damp / smoke is getting out of the pump while the vacuum pump is running, or when the oil is leaked significantly from the element, it is time to change the filter element.

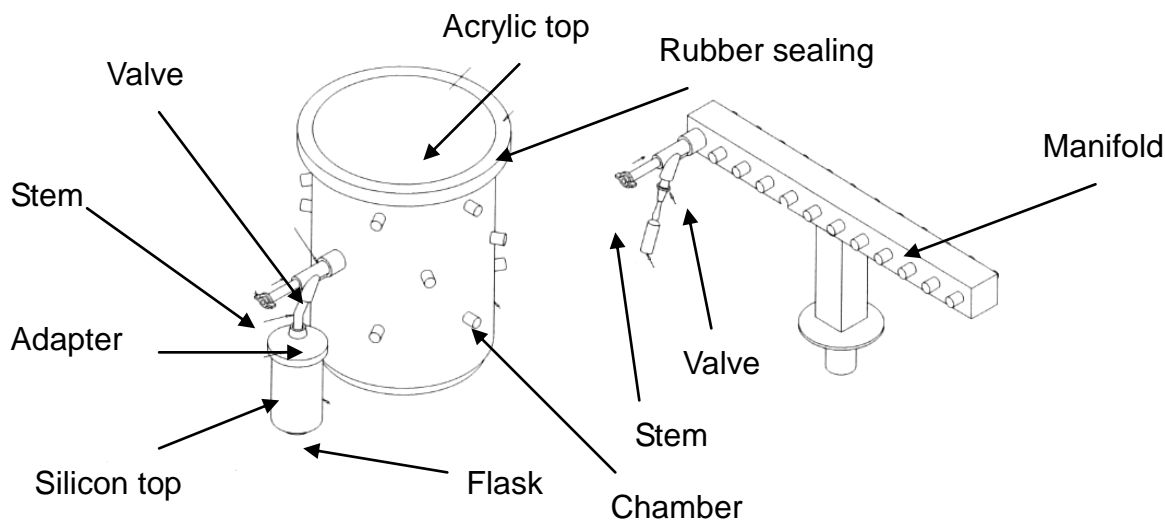
* Remarks: Even though you have recently changed the element, if the smoke and leakage of oil continues, it might be the case that there is a leakage in the vacuum line or the ballast valve is opened too widely. If there is a leakage, please make sure this will be blocked. If the ballast valve is open, please closed it until the smoke disappears.

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6.3. Drying Chamber (드로잉 챔버)


6.3.1. Index of the parts of the drying chamber (부분별 명칭)

6.3.2. How to connect accessories (연결 방법)

[How to connect flask button generally]

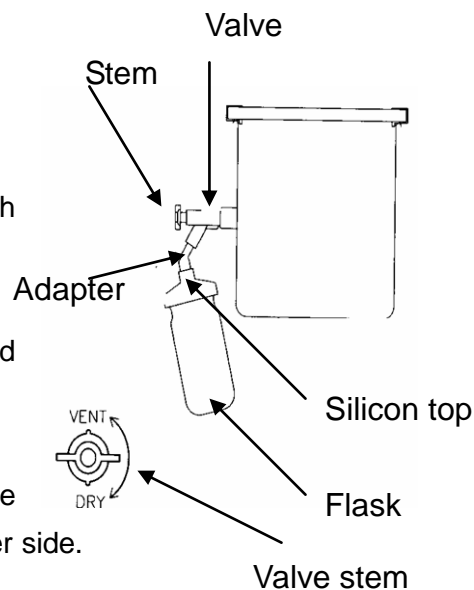- Put the test material on the flask bottom and freeze it covered with plastic wrap, in the Shell Freezer or Deep Freezer under the -30°C .
- Take the plastic wrap of the frozen flask and cover it with the silicon top.
(Please refer the picture alongside)
- Connect the upper part of the silicon top to the adaptor and connect the adaptor to the valve.
(Refer the picture alongside)
- Turn the VENT indicated at the top of the connected valve stem by 180° and adjust that the DRY comes to the upper side.
(Please refer the picture alongside)
- The test sample starts to dry.

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[How to connect ampules generally to the system]

- Put the test sample in the ampule and connect the opening to the ampule adaptor.
- Turn the VENT indicated at the top of the stem by 180° and adjust that the DRY comes to the upper side.
- The test sample starts to dry.

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7. General Operation of the Freeze Dryer

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1. Before you operate the vacuum pump, turn on the freeze dryer first.
2. Please wait until the temperature of the cold trap goes down to -40°C
3. Before you operate the vacuum pump, please check the oil quality and level. If the oil quality is bad, please change the oil.
 - * When the vacuum pressure is under 50mTorr, connect the sample to the valve.
4. To prevent moisturizing in the pump, please, open the gas ballast of the vacuum pump.
5. In the case the operation of the freeze dryer is stopped, after the test:
 - Turn off the vacuum pump
 - Open the valve inside of the chamber to release vacuum (or open the drain valve)
 - Turn the freeze dryer off
 - Defrost the ice layer on the cold trap
 - Expel the moisture from the chamber and wipe out the inside of the chamber, and keep the freezer in clean condition.

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1. When the room temperature is above + 30°C. When the room Temperature is above +30°C, the condenser cannot be cooled. It can be one of cause of trouble.
=> If the room temperature around the unit is +28°C, please often the condenser door and window to optimize the ventilation and cooling.

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2. When the room temperature is under 0°C. The oil of the **compressor/ vacuum pump** is condensed. It can have negative effect on the operation of the **compressor/ vacuum pump**. Each electrical part can be wrongly operated
=> Under the circumstance that the room temperature is under 0°C, make the place warm by using the other heating application and keep the temperature minimal +5°C.

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3. In case there is excessive dust around the unit. This unit use the condenser which is air cooled, In case of the excessive dust around the unit, dust will pill up the front of the condenser. This will cause the compressor to nor work efficiently, This is one of the cause of a defect compressor

=> In case there is excessive dust around the unit door which covers condenser should be often once in a month and remove dust and clean up. Also wipe the dust off the condenser with a soft brush.

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4. Under heavy humid circumstances. It can cause a technical malfunction in the
electrical application and short circuit.

=> Ventilation the place properly by using the ventilator and opening the windows.
An application, which can reduce moisture, will be helpful.

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8.2. Registration of Service (서비스 등록)

* Before you register for the service, please read the following information:

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1. The unit does not operate at all.

○ Check the power supply (230V)

○ Please, check the power supply plug connection.

○ Please, check whether you turn on the unit according the instructions.

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2. The freeze dryer does not work

○ Do not place heat-emitting appliances around the freeze dryer

○ Please, check whether the condenser is blocked by dust.

○ Please, check whether too much test sample is used at one time.

○ Please, check whether the pre-freezing time of the test sample is long enough.

○ Please, check whether you turn on the unit according the instructions.

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3. No vacuum condition

- Please, check whether the vacuum pump works.
- Please, check whether the oil change period of the vacuum pump has passed.
- Please, check whether the drain valve is closed.
- Please, check whether the vacuum valve is closed.
- Please, check whether too much test sample is used at one time.
- Please, check whether the pre-freezing time of the test sample is long enough.
- Please, check whether you turn on the unit according the instructions.

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4. Superfluous oil at the bottom of the unit

○ During operation of the vacuum pump, some oil can flow out of the air outlet of the vacuum pump.

(Remedy: change the element of the exhaust filter at the upper part of the vacuum pump)

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5. Loud running noise

- Please, check whether the floor is even.
- Please, check whether the noise comes from the compressor.
- * The noise of the compressor is bigger than that of household refrigerators.

O Please, check whether the noise comes from the vacuum pump.

* When the noise is bigger than normal, please check the following three things.

- Please, check whether the vacuum pump is installed on the right position.

=>Remedy: in the case of a wrong position, please move the pump a little bit.

- Please, check the vacuum circulation, a valve or the drain valve might not be closed.

=>Remedy: in the case that they are open, please close them.

- Please, check whether the gas ballast located at the upper part of the vacuum pump is open too widely.

=>When the ballast valve is open too widely, please turn it to the right (clockwise) a little bit.

O Do you hear water stream?

* The unit is equipped with many circulation of refrigerant. You can hear the freezing material stream through the circulation. It is not a technical problem.

O Do you hear a cracking sound when you turn on the unit?

* Because the electrical operation of this unit is run by the magnetic relay system, you can hear the cracking sound. However, it is not a technical problem.

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6. Do you smell something unusual?

○ If you smell this while the vacuum pump is running, please, check the following three things

- It can be the case that the residue of old test sample is mixed with the vacuum pump oil.

=> Remedy: Please, change the oil of the vacuum pump.

- It can be the case that the ballast valve is open too widely, and there is a leakage in the

vacuum line.

=>Remedy: Please, close the ballast valve by turning to the right (clockwise) and make the

leakage in the vacuum line closed.

- It can be the case that the element of the exhaust filter should be changed.

=>Please change the element of the exhaust filter.

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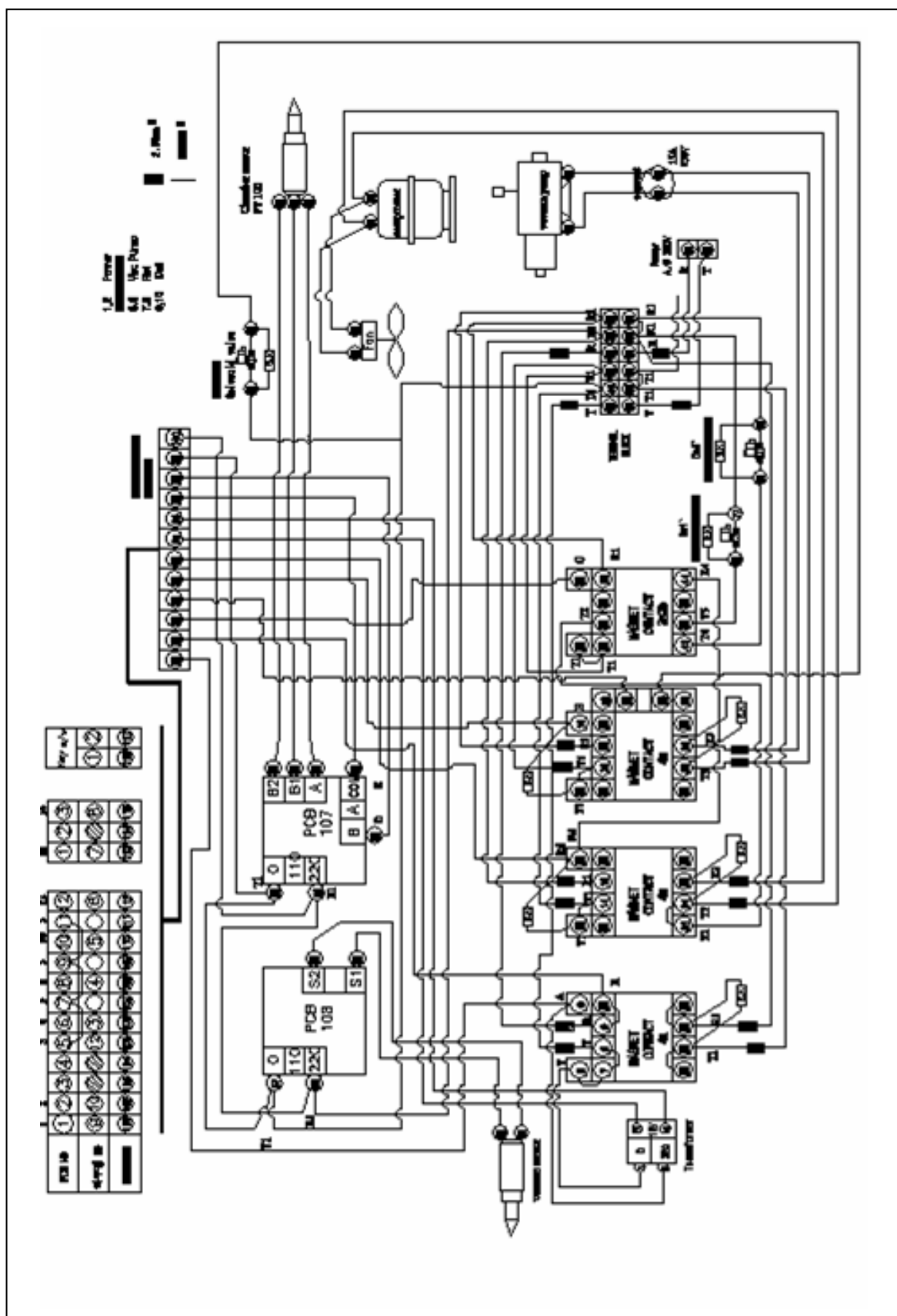
Warranty and contact information

International Warranty

The warranty period starts on the moment the unit leaves the production facility. After this date one year warranty on the parts is valid. During this first year component parts proven to be defective could be send to your local dealer and will be covered by the first year warranty on parts only. Expendable items, i.e., glass, filters, pilot lights, light bulbs and door gaskets are excluded from this warranty. IIShinbiobase Co., Ltd shall not be liable for any indirect or consequential damages including, without limitation, damages relating to lost profits or loss of products.

Conditions and terms of sale

All prices are ex works from the IIShinbiobase Co. Ltd Factory in Korea, excluding insurance, seaworthy packing and customs duty. Effective prices are those valid on date of confirmation of order. All specifications mentioned in the price list are based on 230 V, 50/60Hz power supply. Continuing research and improvements may result in specification changes. Errors and misprints excepted

Contact information

In case service of sales back up is needed please do not hesitate to contact your local IIShin Dealer:

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Or contact the IIShinbiobase Co., Ltd Headquarter: